

Installation manual Coffalyser.Net

| | |
|---|-----------|
| Introduction | 2 |
| Prerequisites | 2 |
| Default installation - SQL server not present | 3 |
| Procedure default installation Coffalyser.Net | 3 |
| Installation Coffalyser.Net - SQL server present | 15 |
| Procedure installation Coffalyser.Net - SQL server present | 15 |
| License | 23 |
| Procedure install license by downloading from within Coffalyser.Net | 23 |
| Procedure install license by importing license file | 25 |
| Updating Coffalyser.Net | 27 |
| Procedure uninstalling a previous version of Coffalyser.Net | 28 |
| Procedure installing new version – preconfigured SQL server | 30 |
| Procedure installing new version – not preconfigured SQL server | 37 |
| More information & contact | 44 |

Introduction

Coffalyser.Net consists of 3 components that interact with each other:

1. The Coffalyser.Net Client. This is the end-user application and forms the user interface.
2. The Coffalyser.Net database. All information is stored in the database, including raw data, experiments and user accounts. The database can run on Microsoft SQL Server 2008 and higher and on any computer in a network.
3. The Coffalyser.Net Server application. This application can communicate with both the Coffalyser.Net Client and the Coffalyser.Net database, thereby connecting these components with each other.

For the most common setups, the installation procedure of Coffalyser.Net is straight forward and does not require advanced IT knowledge. The software can be installed on 32 and 64 bit Windows 7, 8 and 10 systems.

For most users the default installation method will suffice (see **Default installation - SQL server not present** from page 3 onwards). However, Coffalyser.Net allows the use of an already present SQL database server (see **Installation Coffalyser.Net - SQL server present** from page 15 onwards).

In case you want to update Coffalyser.Net to the latest available version, please follow the procedures described in the chapter **Updating Coffalyser.Net** (from page 27 onwards)

Prerequisites

For a successful installation of Coffalyser.Net, the following requirements have to be met:

- A genuine copy of Windows is installed
- The latest Windows updates are downloaded and installed
- Microsoft.Net Framework 3.5 is installed
- Windows Installer 4.5 or higher is installed
- 512 MB of RAM is available
- 1 GB free disk space is available
- You have administrative rights (not required for running Coffalyser.Net)

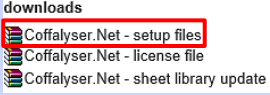
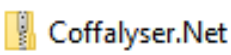
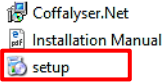
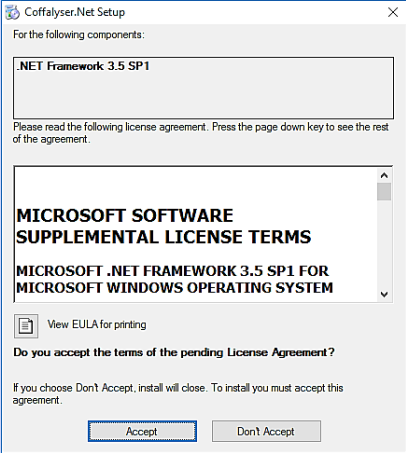
Antivirus software and firewalls might interfere with the installation process. In case the installation of Coffalyser.Net cannot be completed, we recommend to switch these off temporarily and to retry the installation of Coffalyser.Net.

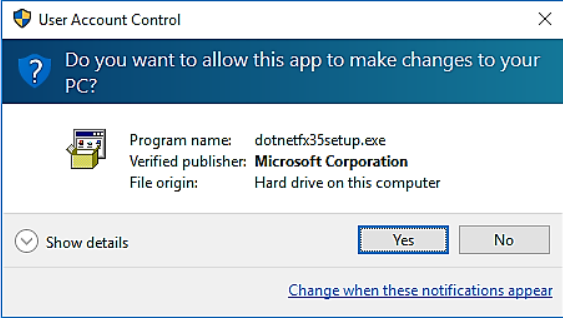
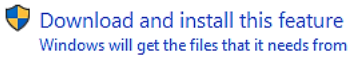
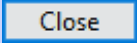
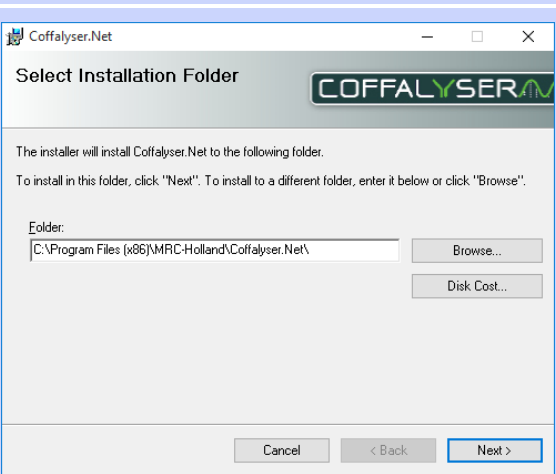
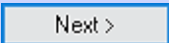


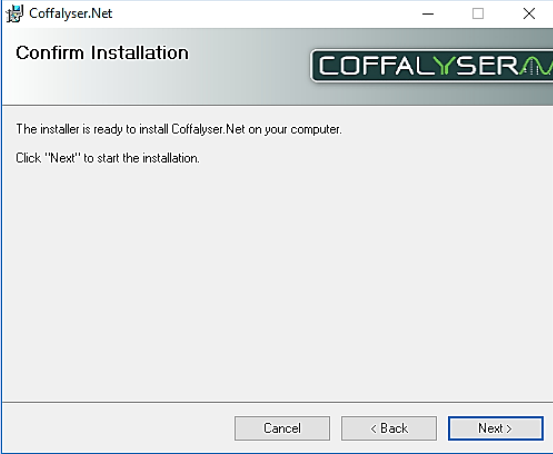
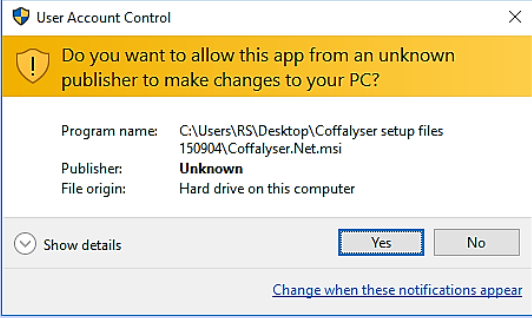
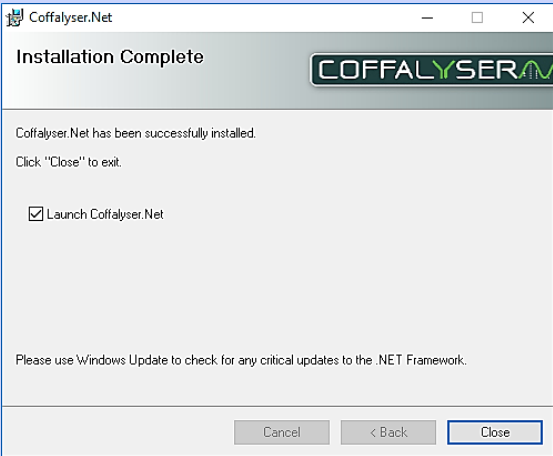
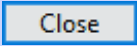
Default installation - SQL server not present


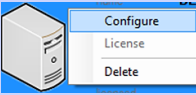
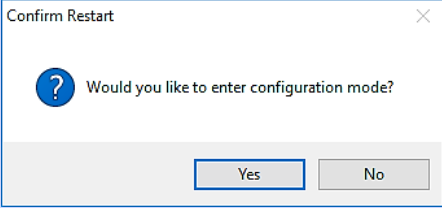
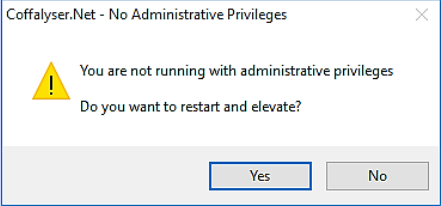
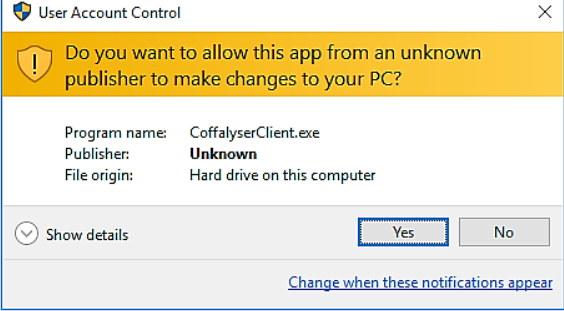
The procedure in this section describes the default installation method for computers on which Coffalyser.Net has not been installed before. This is the recommended method for most users. This method will install all three components (the Client, the database and the Server application) onto the same computer. In case you want to update Coffalyser.Net, please follow the procedures described in the chapter **Updating Coffalyser.Net** from page 27 onwards.

PROCEDURE DEFAULT INSTALLATION COFFALYSER.NET

| | |
|---|---|
| 1. Download the zip file with the Coffalyser.Net – setup files from your MyMLPA account on www.mlpa.com |  |
| 2. Extract the zip file |  |
| 3. Start the installation by opening the setup file |  |
| When prompted with the license agreement for .NET Framework 3.5 SP1, follow steps 4 to 7. Otherwise proceed with step 8. | |
| 4.  Click Accept to accept the license agreement | |

| | | |
|---|---|---|
| 5. |  <p>Program name: dotnetfx35setup.exe Verified publisher: Microsoft Corporation File origin: Hard drive on this computer</p> <p>Yes No</p> <p>Change when these notifications appear</p> | |
| If prompted, click Yes in the User Account Control window | | |
| 6. | Click Download and install this feature |  |
| The download and installation procedure starts. This may take a while. | | |
| 7. | Click Close to finish the installation of .NET Framework 3.5 and continue with the installation of Coffalyser.Net |  |
| The Coffalyser.Net installation wizard opens | | |
| 8. | Maximize the Coffalyser.Net installation wizard from the task bar if this doesn't happen automatically | |
| 9. |  <p>The installer will install Coffalyser.Net to the following folder. To install in this folder, click "Next". To install to a different folder, enter it below or click "Browse".</p> <p>Folder: <input type="text" value="C:\Program Files (x86)\MRC-Holland\Coffalyser.Net\"/> Browse... Disk Cost...</p> <p>Cancel < Back Next ></p> | |
| If it is desired to install Coffalyser.Net on another location than the default, click Browse . If the default location is ok, continue with the next step | | |
| 10. | Click Next |  |

| | | |
|-----------------------|---|---|
| 11. |  | |
| 12. | <p>Click Next to start the installation</p>  <p>If prompted, click Yes in the User Account Control window</p> | |
| 13. |  <p>Ensure that the selection box <i>Launch Coffalyser.Net</i> is checked</p> | |
| 14. | <p>Click Close to exit the installation window</p> |  |
| Coffalyser.Net starts | | |

| | | |
|--|--|---|
| 15. |  <p>License Agreement</p> <p>Coffalyser</p> <p>License</p> <p>Coffalyser End User License Agreement</p> <p>NOTICE TO USER: PLEASE READ THIS CONTRACT CAREFULLY. BY USING ALL OR ANY PORTION OF THE SOFTWARE YOU ACCEPT ALL THE TERMS AND CONDITIONS OF THIS AGREEMENT. YOU AGREE THAT THIS AGREEMENT IS ENFORCEABLE LIKE ANY WRITTEN NEGOTIATED AGREEMENT SIGNED BY YOU. IF YOU DO NOT AGREE, DO NOT USE THIS SOFTWARE. IF YOU ACQUIRED THE SOFTWARE ON TANGIBLE MEDIA (e.g. CD) WITHOUT AN OPPORTUNITY TO REVIEW THIS LICENSE AND YOU DO NOT ACCEPT THIS AGREEMENT: (A) DO NOT USE THE SOFTWARE AND (B) RETURN IT.</p> <p>1. Definitions: "Software" means (a) all of the contents of the files, disk(s), CD-ROM(s) or other media with which this Agreement is provided, including but not limited to (i) Coffalyser (MRC-Holland), or third party computer information or software; (ii) digital images, stock photographs, clip art, sounds or other artistic works ("Stock Files"); (iii) related explanatory written materials or files ("Documentation"); and (iv) fonts; and (b) upgrades, modified versions, updates.</p> <p>Please read the agreement above carefully and choose one of the options below</p> <p><input type="button" value="I Accept"/> <input type="button" value="I Do Not Accept"/></p> | |
| 16. | <p>In the Coffalyser.Net Server Selection window right mouse click on the (grey) server with the name of your computer and select <i>Configure</i></p> |  |
| 17. |  <p>Confirm Restart</p> <p>Would you like to enter configuration mode?</p> <p><input type="button" value="Yes"/> <input type="button" value="No"/></p> | |
| 18. |  <p>Coffalyser.Net - No Administrative Privileges</p> <p>You are not running with administrative privileges Do you want to restart and elevate?</p> <p><input type="button" value="Yes"/> <input type="button" value="No"/></p> | |
| 19. |  <p>User Account Control</p> <p>Do you want to allow this app from an unknown publisher to make changes to your PC?</p> <p>Program name: CoffalyserClient.exe Publisher: Unknown File origin: Hard drive on this computer</p> <p><input type="button" value="Yes"/> <input type="button" value="No"/></p> <p>Change when these notifications appear</p> | |
| <p>The Coffalyser.Net – Configuration Wizard opens</p> | | |

20.



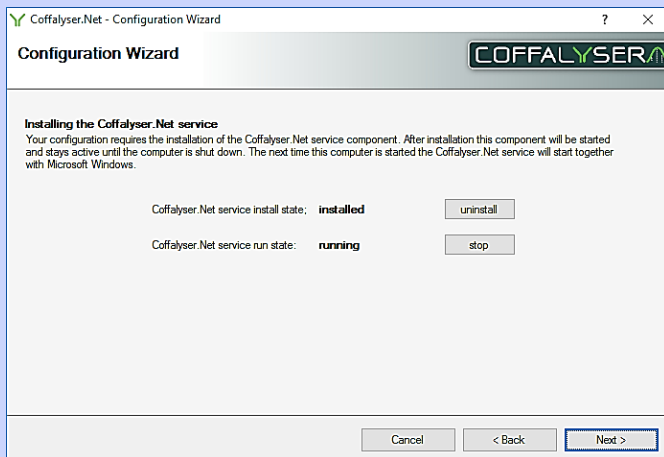
Click **Next**

21.




Click **install** to install the Coffalyser.Net service

22.





Click **Next**

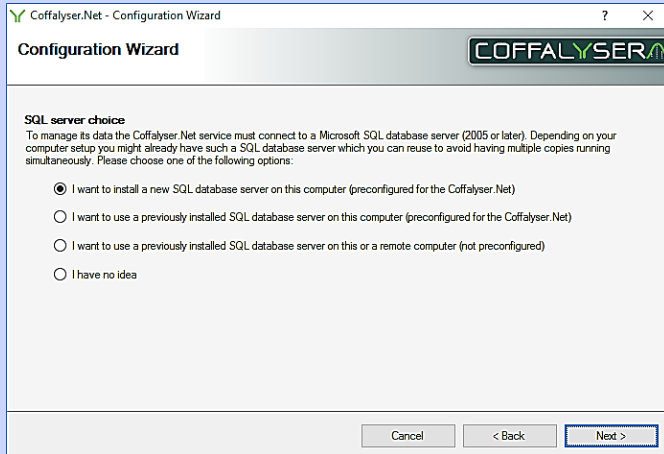

23.

Select the option *I want to install a new SQL database server on this computer (preconfigured for the Coffalyser.Net)*

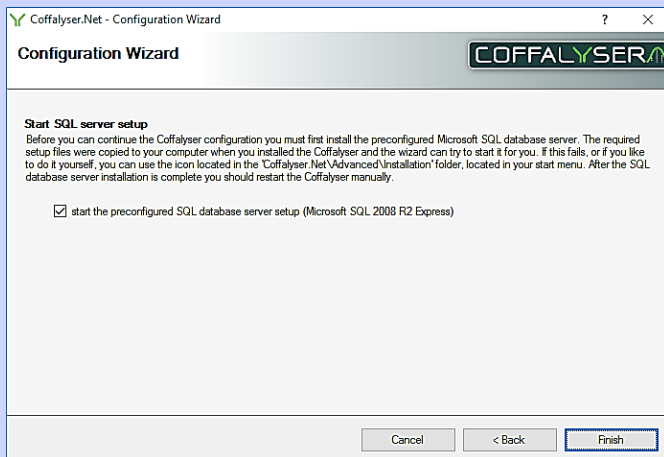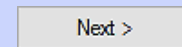
24. Click **Next**



25.

Ensure that the selection box *start the preconfigured SQL database server setup (Microsoft SQL 2008 R2 Express)* is checked

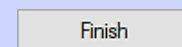26. Click **Finish**



The SQL Server 2008 R2 Express installation wizard opens

27. Maximize the SQL Server 2008 R2 Express installation wizard from the task bar if this doesn't happen automatically

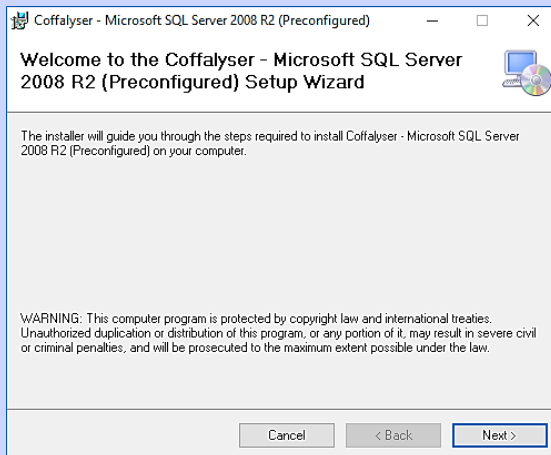
28.



Click **Accept** to accept the license agreement


The download and installation procedure starts. This may take a while.

29.






Click **Next**

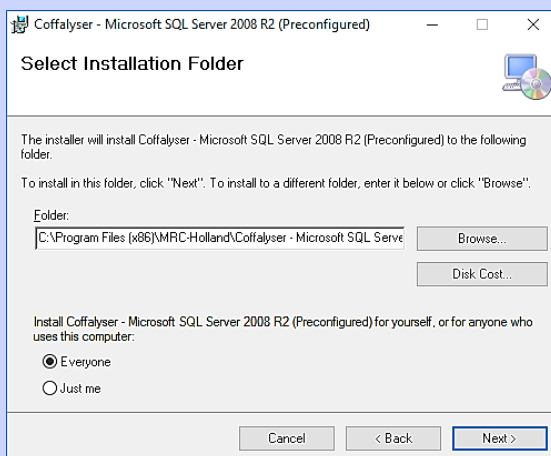
30.



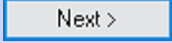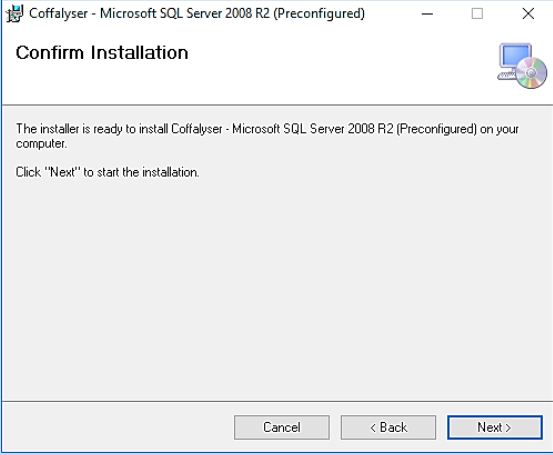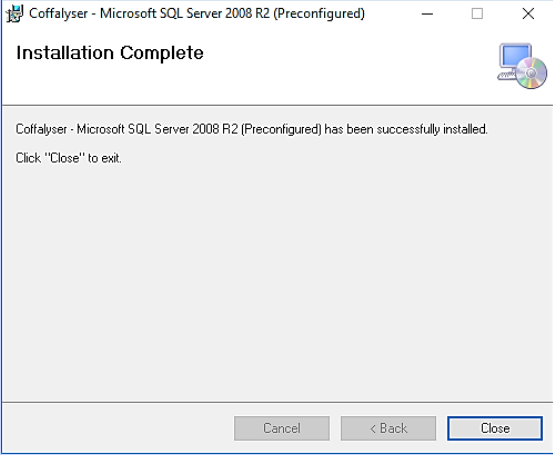Ensure that the option *Everyone* is selected


| | |
|---|--|
| 31. Click Next |  |
| 32. |  <p>The installer is ready to install Coffalyser - Microsoft SQL Server 2008 R2 (Preconfigured) on your computer. Click "Next" to start the installation.</p> <p>Buttons: Cancel, < Back, Next ></p> <p>Click Next to start the installation</p> |
| 33. |  <p>Coffalyser - Microsoft SQL Server 2008 R2 (Preconfigured) has been successfully installed. Click "Close" to exit.</p> <p>Buttons: Cancel, < Back, Close</p> <p>Click Close to exit the installation wizard</p> |
| 34. Double click on the Coffalyser.Net shortcut icon on the desktop to start the software |  |
| 35. In the Coffalyser.Net Server Selection window right mouse click on the (orange) server with the name of your computer and select <i>Configure</i> |  |
| 36. |  <p>Would you like to enter configuration mode?</p> <p>Buttons: Yes, No</p> <p>Click Yes to enter the configuration mode</p> |

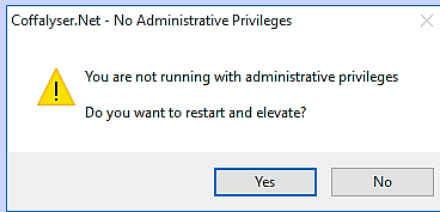37.



If prompted, click **Yes** to restart and elevate in order to run with administrative privileges

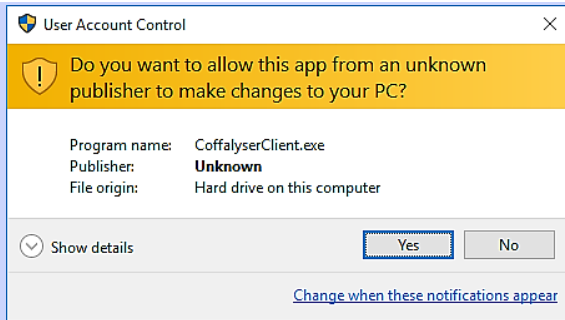
38.



If prompted, click **Yes** in the User Account Control window


The Coffalyser.Net – Configuration Wizard opens

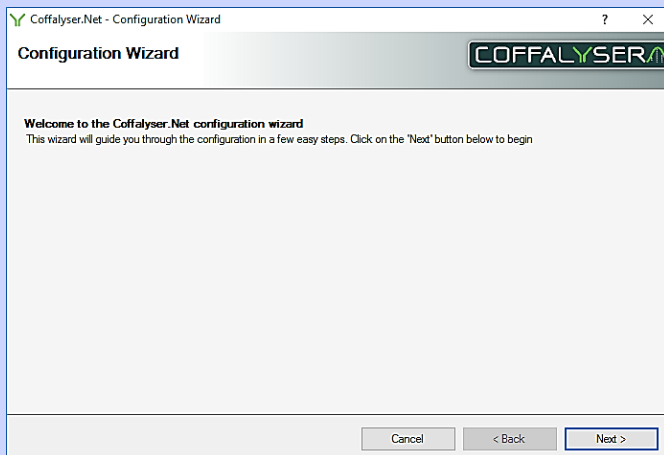39.




Click **Next**

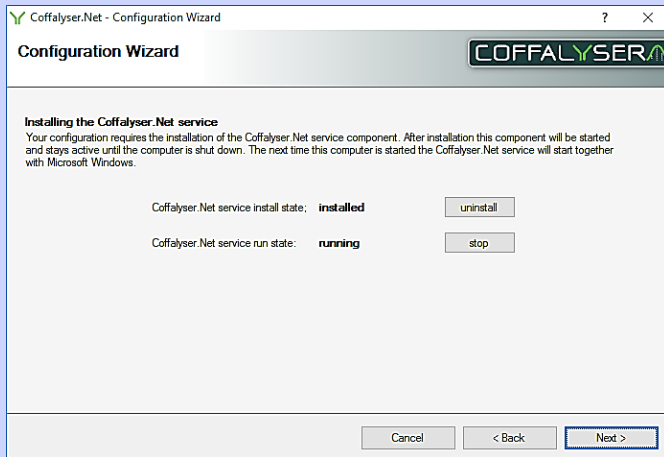40.




Click **Next**

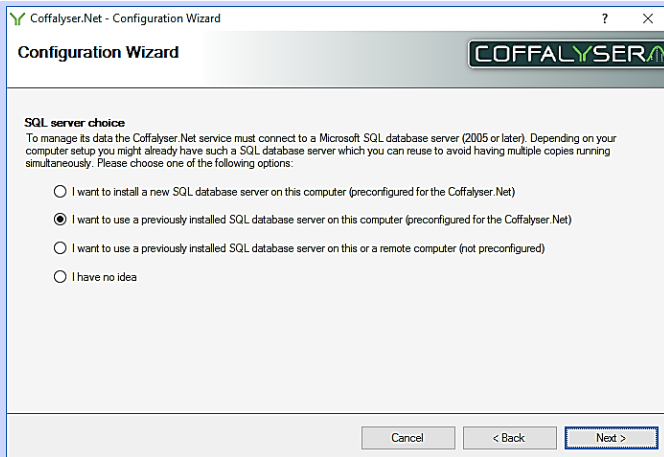41.




Select the option *I want to use a previously installed SQL database server on this computer (preconfigured for the Coffalyser.Net)* if it is not preselected

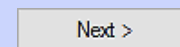42. Click **Next**



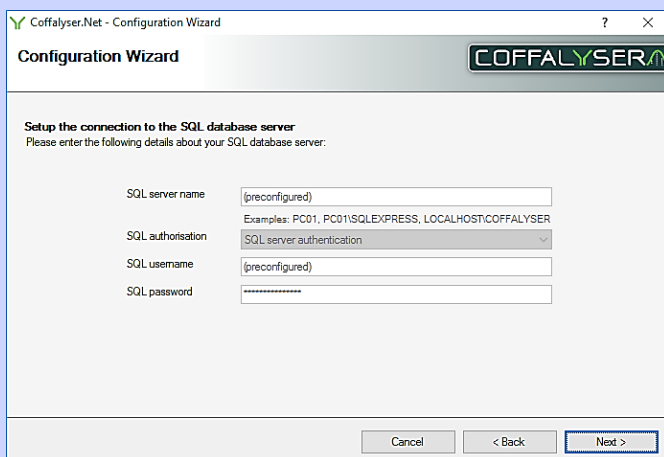43.



Click **Next**

44.

When desired, change the name of the database in the designated field. By default it is Coffalyser

45. Click **Next**

Next >

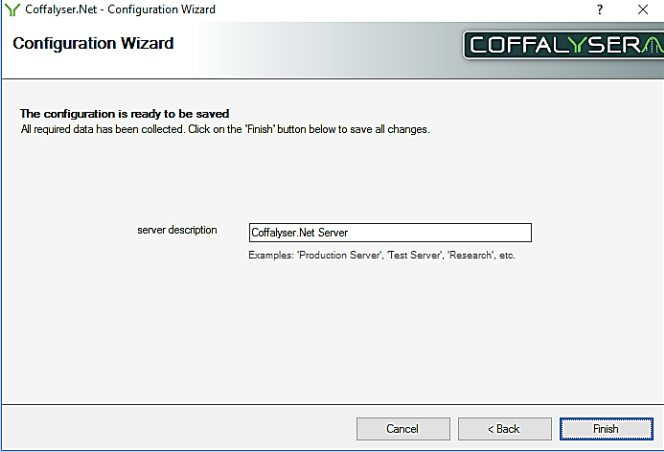
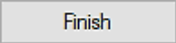
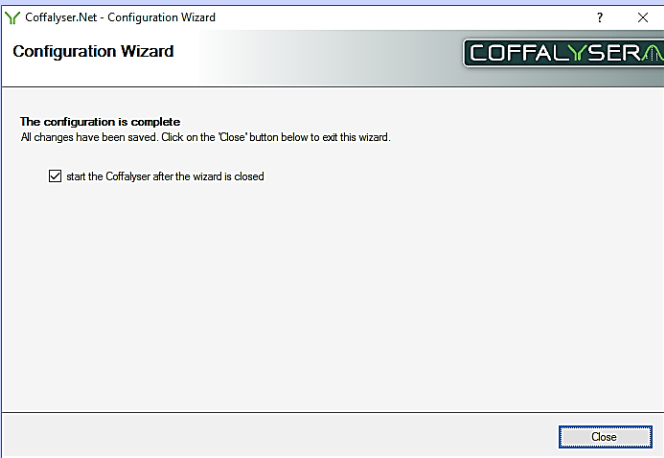
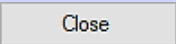
46.

Create a name for the first organisation, a password, a secret question and answer. Enter these in the respective fields

Write down or remember the entered password, secret question and answer. These are needed later to log in to Coffalyser.Net.

47. Click **Next**





Next >






| | | |
|-----|--|---|
| 48. |  <p>When desired, change the name of the server in the designated field. By default it is Coffalyser.Net Server</p> | |
| 49. | Click Finish |  |
| 50. |  <p>Ensure that the selection box <i>start the Coffalyser after the wizard is closed</i> is checked</p> | |
| 51. | Click Close to exit the Configuration wizard |  |
| 52. | Install your license as described in section in the chapter License on page 23 | |


Installation Coffalyser.Net - SQL server present

The procedure in this section describes the installation method for Coffalyser.Net when an SQL server is already present on the computer and on which Coffalyser.Net has not been installed before. Please note that this method requires specific details of the SQL server (e.g. location, login credentials, authentication method etc.). If you do not know these, it is recommended to use the default installation method for Coffalyser.Net (see the previous section). In case you want to update or reinstall Coffalyser.Net, please follow the procedures described in the chapter **Updating Coffalyser.Net** from page 27 onwards

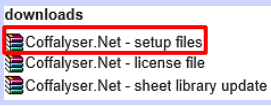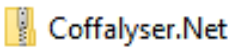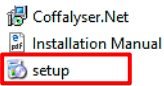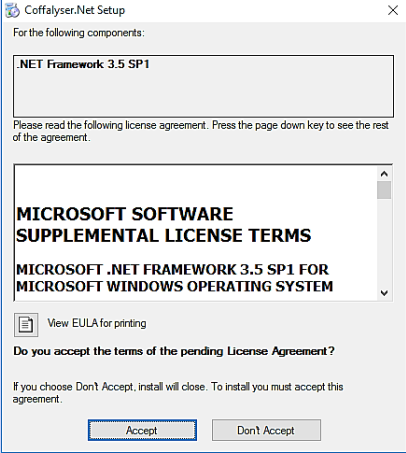
PROCEDURE INSTALLATION COFFALYSER.NET - SQL SERVER PRESENT

| | |
|--|---|
| <p>1. Download the zip file with the Coffalyser.Net – setup files from your MyMLPA account on www.mlpa.com</p> |  |
| <p>2. Extract the zip file</p> |  |
| <p>3. Start the installation by opening the setup file.</p> |  |
| <p>When prompted with the license agreement for .NET Framework 3.5 SP1, follow steps 4 to 7. Otherwise proceed with step 8.</p> | |
| <p>4.</p>  <p>Click Accept to accept the license agreement</p> | |

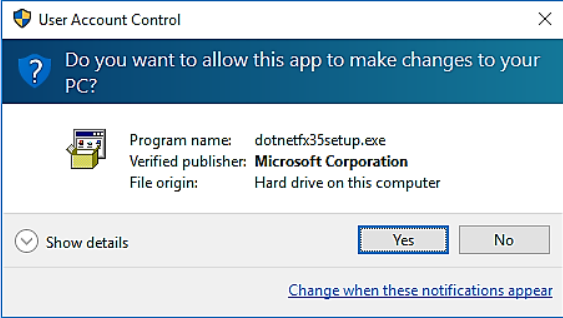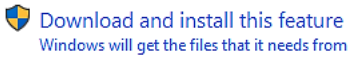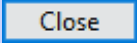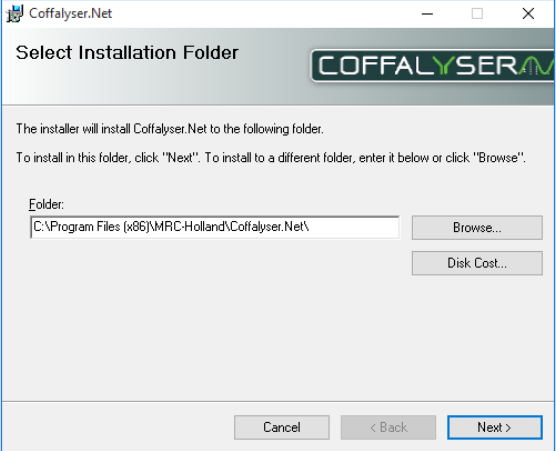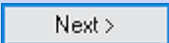
| | | |
|---|---|---|
| 5. |  <p>The dialog box asks: "Do you want to allow this app to make changes to your PC?". It shows details for the program: "Program name: dotnetfx35setup.exe", "Verified publisher: Microsoft Corporation", and "File origin: Hard drive on this computer". There are "Yes" and "No" buttons, and a "Show details" link.</p> | |
| If prompted, click Yes in the User Account Control window | | |
| 6. | Click Download and install this feature |  |
| The download and installation procedure starts. This may take a while. | | |
| 7. | Click Close to finish the installation of .NET Framework 3.5 and continue with the installation of Coffalyser.Net |  |
| The Coffalyser.Net installation wizard opens | | |
| 8. | Maximize the Coffalyser.Net installation wizard from the task bar if this doesn't happen automatically. | |
| 9. |  <p>The dialog shows the default folder: "C:\Program Files (x86)\MRC-Holland\Coffalyser.Net\". There are "Browse...", "Disk Cost...", "Cancel", "< Back", and "Next >" buttons.</p> | |
| If it is desired to install Coffalyser.Net on another location than the default, click Browse . If the default location is ok, continue with the next step | | |
| 10. | Click Next |  |

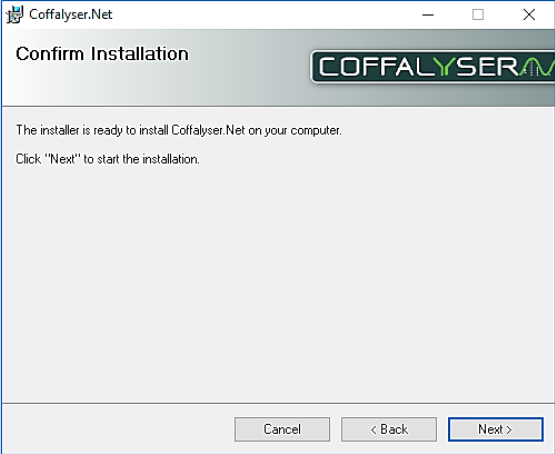11. 

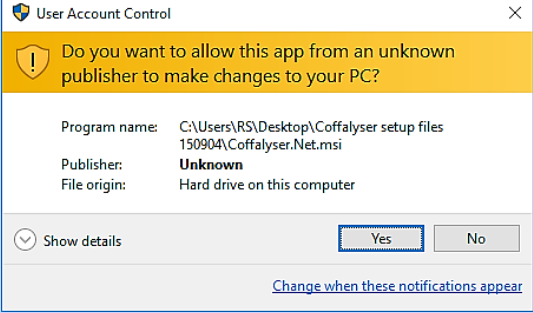Click **Next** to start the installation

12. 

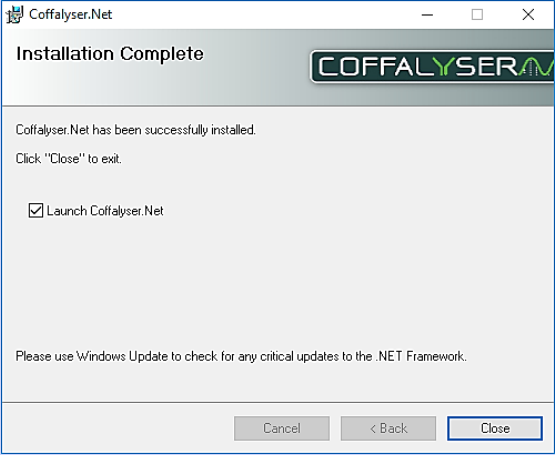If prompted, click **Yes** in the User Account Control window

13. 

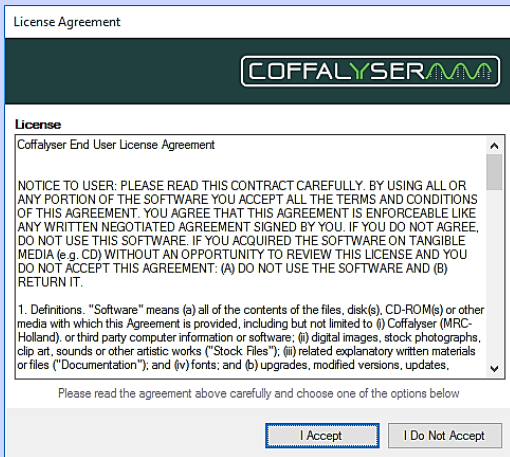
Ensure that the selection box *Launch Coffalyser.Net* is checked

14. Click **Close** to exit the installation window



Coffalyser.Net starts







15.


Click **I Accept** to accept the license agreement of Coffalyser.Net

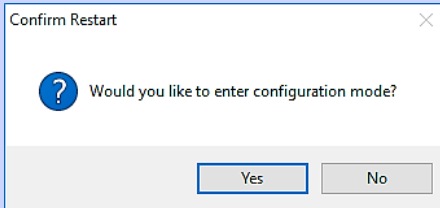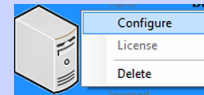16. In the Coffalyser.Net Server Selection window right mouse click on the (grey) server with the name of your computer and select *Configure*




17.

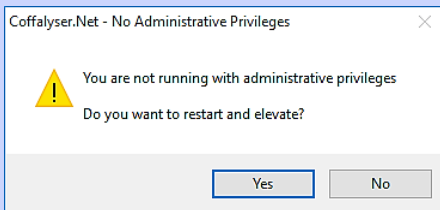Click **Yes** to enter the configuration mode



18.

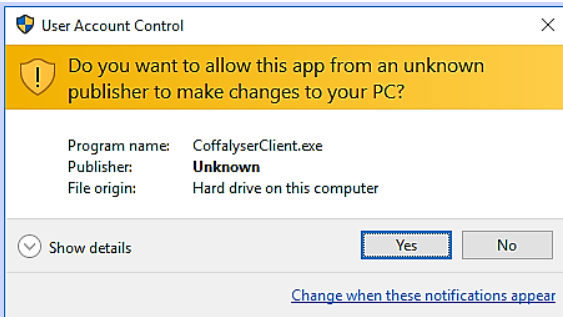If prompted, click **Yes** to restart and elevate in order to run with administrative privileges




19.

If prompted, click **Yes** in the User Account Control window


The Coffalyser.Net – Configuration Wizard opens

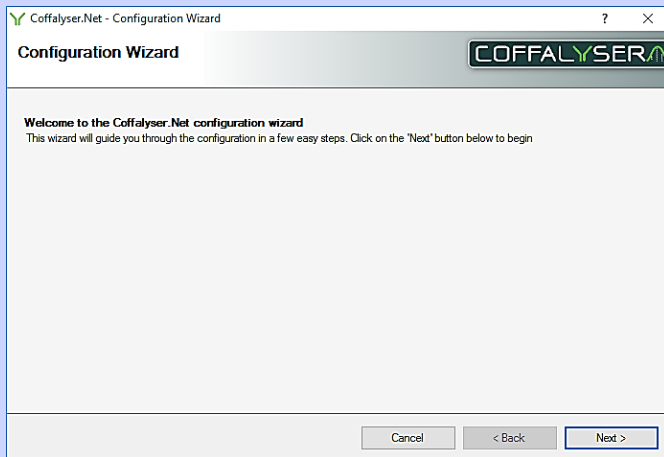20.




Click **Next**

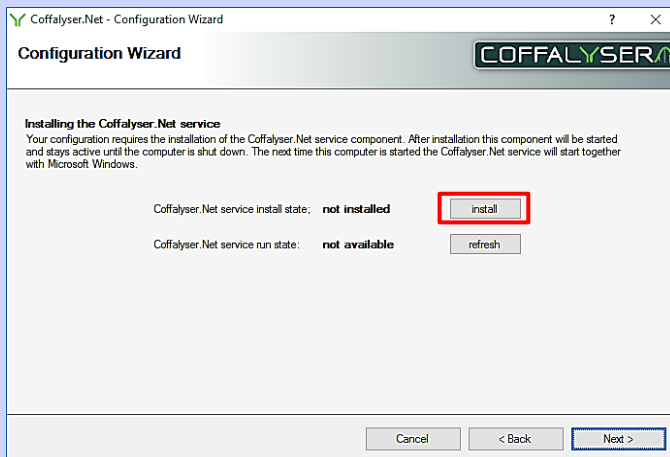21.




Click **install** to install the Coffalyser.Net service

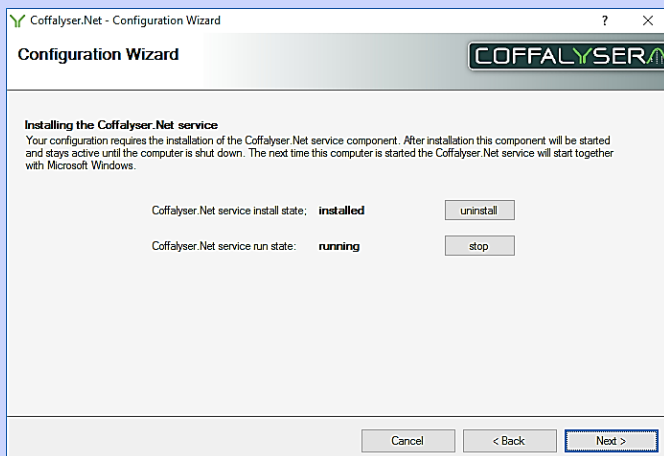22.





Click **Next**

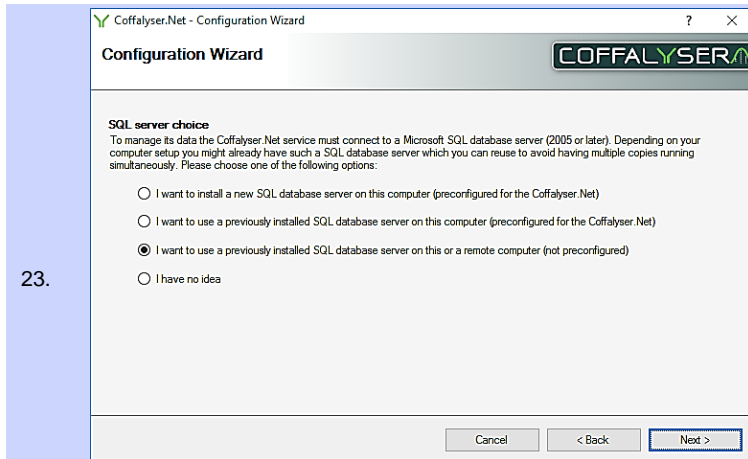


Select the option *I want to use a previously installed SQL database server on this or a remote computer (not preconfigured)*

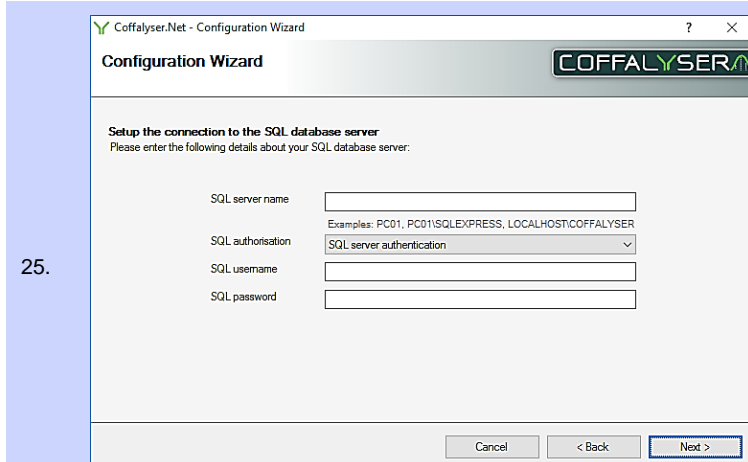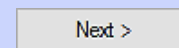24. Click **Next**



Enter the information of the SQL server in the designated fields

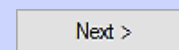
26. Click **Next**



27.

When desired, change the name for the database in the designated field. By default it is Coffalyser

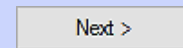28. Click **Next**








29.

Create a name for the first organisation, a password, a secret question and answer. Enter these in the respective fields

Write down or remember the entered password, secret question and answer. These are needed later to log into Coffalyser.Net.

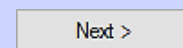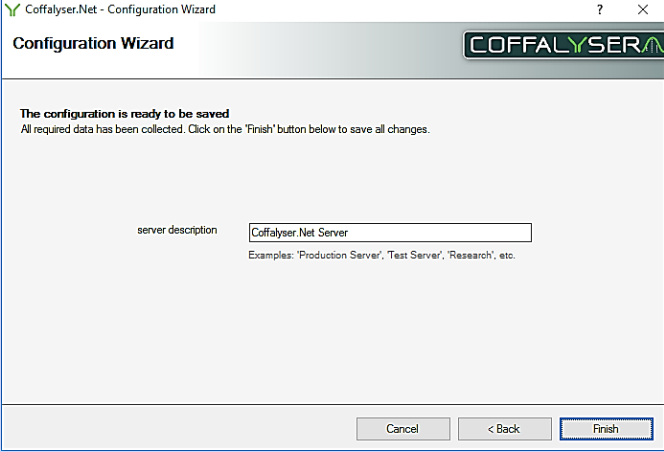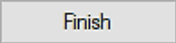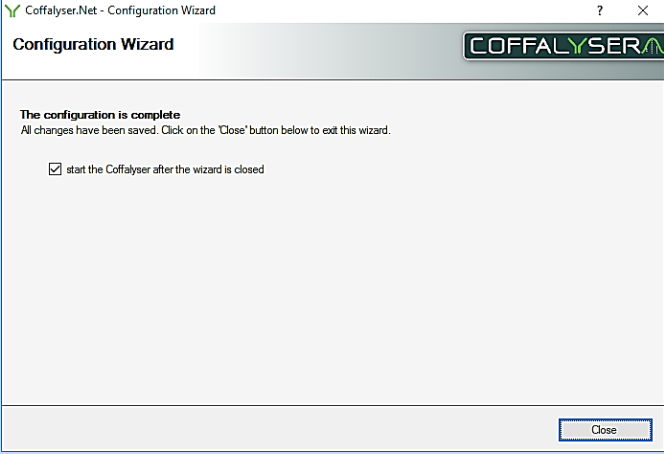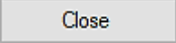30. Click **Next**


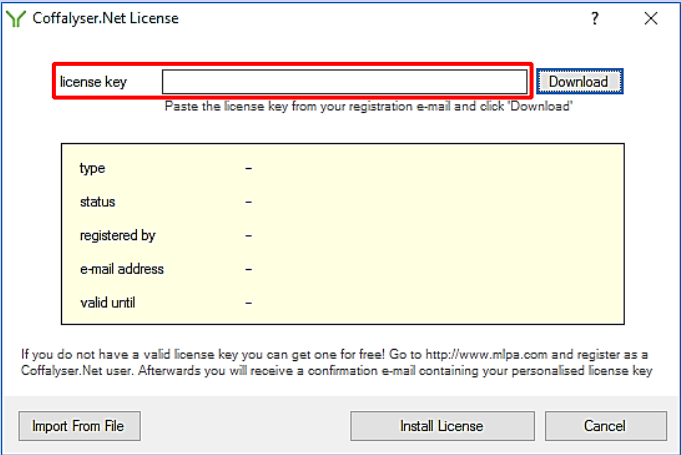
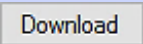
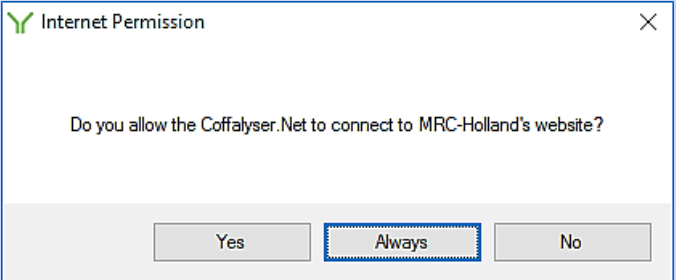


| | | |
|-----|--|---|
| 31. |  <p>When desired, change the name of the server in the designated field. By default it is Coffalyser.Net Server</p> | |
| 32. | Click Finish |  |
| 33. |  <p>Ensure that the selection box <i>start the Coffalyser after the wizard is closed</i> is checked</p> | |
| 34. | Click Close to exit the Configuration wizard |  |
| 35. | Install your license as described in section in the chapter License on page 23 | |


License

Before you can log in and use Coffalyser.Net, it is necessary to activate the software by installing your license key. This can be done by downloading the license directly from within Coffalyser.Net. When you cannot download the license (because you do not have internet access, for instance), you can also import the license file. This file can be downloaded from your MyMLPA account on www.mlpa.com.

PROCEDURE INSTALL LICENSE BY DOWNLOADING FROM WITHIN COFFALYSER.NET


| | |
|--|---|
| 1. In the Coffalyser.Net Server Selection window right mouse click on the (orange) server with the name of your computer and select <i>License</i> |  |
| 2.  Enter your license key in the designated field. ¹ | |
| 3. Click Download |  |
| 4.  Click Always to allow Coffalyser.Net to connect to the MRC-Holland server to download the license | |

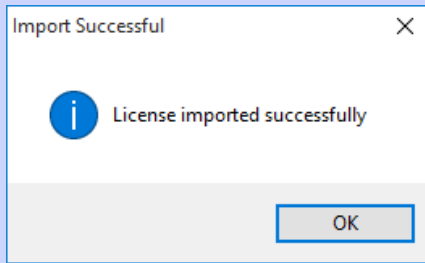¹ Your license key has been sent to you by e-mail. It can also be found in the Coffalyser.Net section in your MyMLPA account on www.mlpa.com.



5.

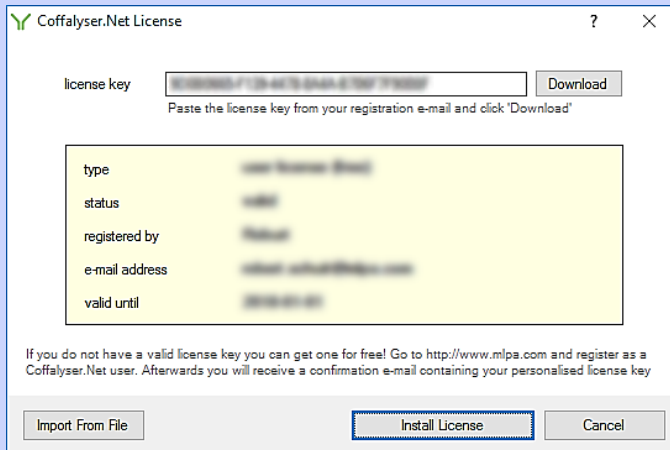Click **OK**











6.

Click **Install License**



7. Coffalyser.Net is now ready for use

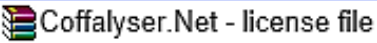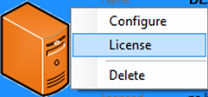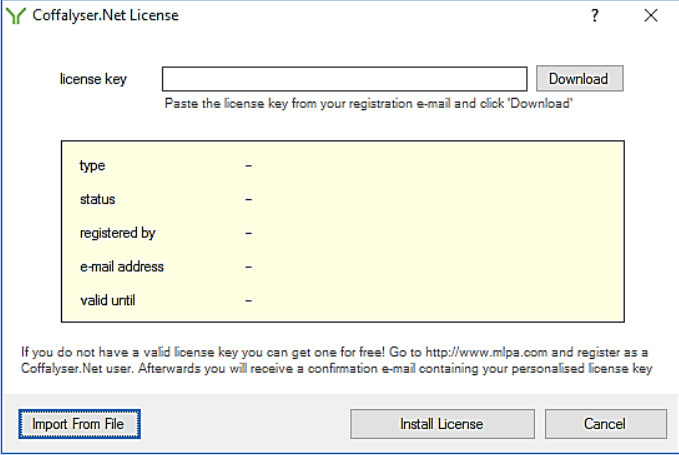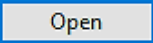PROCEDURE INSTALL LICENSE BY IMPORTING LICENSE FILE

| | |
|---|---|
| <p>1. Log in to your MyMLPA account on www.mlpa.com</p> |  |
| <p>2. Navigate to the section my Coffalyser.Net</p> |  |
| <p>3. Click on Coffalyser.Net – license file to download the Coffalyser.license file</p> |  |
| <p>4. Save the file on your computer</p> | |
| <p>If you downloaded the license file on another computer than on which Coffalyser.Net has been installed, copy this file to an USB drive. Next, insert the USB drive in the computer on which Coffalyser.Net has been installed.</p> | |
| <p>5. In the Coffalyser.Net Server Selection window right mouse click on the (orange) server with the name of your computer and select <i>License</i></p> |  |
| <p>6. </p> <p>Click Import From File</p> | |
| <p>A dialog box opens</p> | |
| <p>7. Navigate to the location of the Coffalyser.license file</p> | |
| <p>8. Select the file and click Open</p> |  |

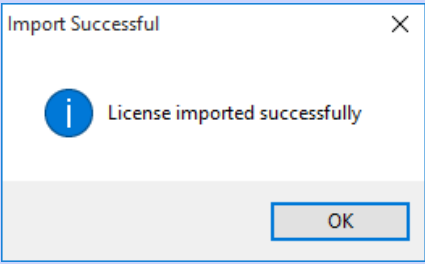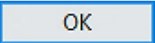9.  

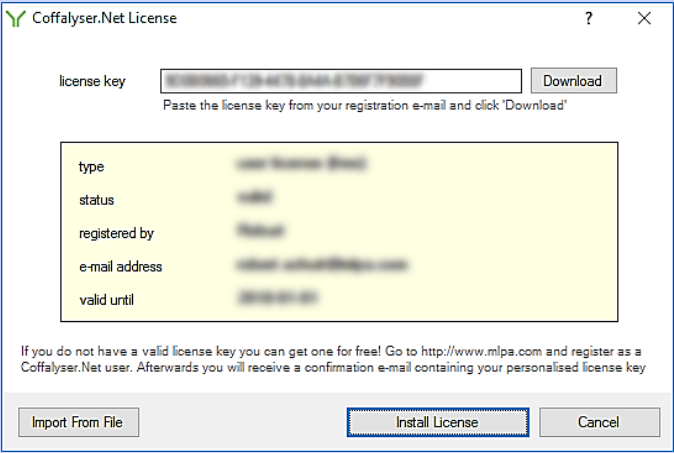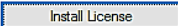Click **OK**

10.  

Click **Install License**

11. Coffalyser.Net is now ready for use

Updating Coffalyser.Net

Based on user feedback, new insights and/or regulations, Coffalyser.Net is continuously improved. This results in new versions of the software, which are regularly released.

Coffalyser.Net can easily be updated to the latest version by uninstalling the previous version and installing the latest one.

Important notes:


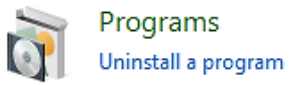
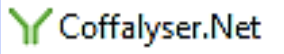
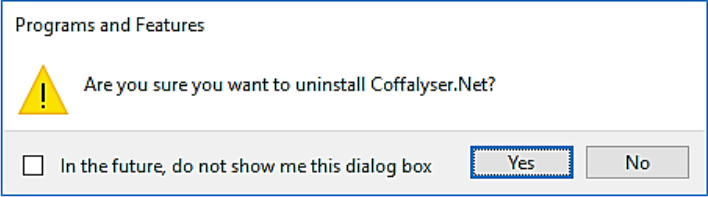
- **Do not uninstall any version of Microsoft SQL server. This also include Coffalyser - Microsoft SQL Server 2008 R2 (preconfigured)!**



As Microsoft SQL Server host the database of Coffalyser.Net, uninstalling this program may lead to irreversible data loss!

- For installation of the new version of Coffalyser.Net it is important that the correct procedure is followed:
 - When for the original installation a preconfigured SQL database server was used, see **Procedure installing new version – preconfigured SQL server**
 - When for the original installation a previously installed, not preconfigured SQL database server was used, see **Procedure installing new version – not preconfigured SQL server**

PROCEDURE UNINSTALLING A PREVIOUS VERSION OF COFFALYSER.NET






This procedure has been described based on the standard configuration in Windows 10. The location of the Control Panel and the option 'Uninstall a program' may be different in other Windows versions and other configurations. If you are unable to find these, consult the documentation on the operating system or contact the IT department of your institute for assistance.





| 1. Open the Control Panel of Windows |  | | | | | | | | | | |
|---|---|-----------|---|-------------|-----------------------|-------------|---------------------------------------|-----------------------|--|-----------------------|---|
| 2. Click Uninstall a program in the category PROGRAMS |  | | | | | | | | | | |
| 3. <table border="1" data-bbox="256 730 984 903"><thead><tr><th>Name</th><th>Publisher</th></tr></thead><tbody><tr><td>Coffalyser - Microsoft SQL Server 2008 R2 (Preconfig...</td><td>MRC-Holland</td></tr><tr><td>Coffalyser.Net</td><td>MRC-Holland</td></tr><tr><td>Microsoft SQL Server 2008 R2 (64-bit)</td><td>Microsoft Corporation</td></tr><tr><td>Microsoft SQL Server 2008 R2 Native Client</td><td>Microsoft Corporation</td></tr></tbody></table> Select <i>Coffalyser.Net</i> from the list | Name | Publisher | Coffalyser - Microsoft SQL Server 2008 R2 (Preconfig... | MRC-Holland | Coffalyser.Net | MRC-Holland | Microsoft SQL Server 2008 R2 (64-bit) | Microsoft Corporation | Microsoft SQL Server 2008 R2 Native Client | Microsoft Corporation |  |
| Name | Publisher | | | | | | | | | | |
| Coffalyser - Microsoft SQL Server 2008 R2 (Preconfig... | MRC-Holland | | | | | | | | | | |
| Coffalyser.Net | MRC-Holland | | | | | | | | | | |
| Microsoft SQL Server 2008 R2 (64-bit) | Microsoft Corporation | | | | | | | | | | |
| Microsoft SQL Server 2008 R2 Native Client | Microsoft Corporation | | | | | | | | | | |
| WARNING: Do <u>NOT</u> select any version of Microsoft SQL Server. Uninstalling may lead to irreversible data loss! | | | | | | | | | | | |
| 4. <table border="1" data-bbox="256 1077 984 1255"><thead><tr><th>Name</th><th>Publisher</th></tr></thead><tbody><tr><td>Coffalyser - Microsoft SQL Server 2008 R2 (Preconfig...</td><td>MRC-Holland</td></tr><tr><td>Coffalyser.Net</td><td>MRC-Holland</td></tr><tr><td>Microsoft SQL Server 2008 R2 (64-bit)</td><td>Microsoft Corporation</td></tr></tbody></table> Click Uninstall | Name | Publisher | Coffalyser - Microsoft SQL Server 2008 R2 (Preconfig... | MRC-Holland | Coffalyser.Net | MRC-Holland | Microsoft SQL Server 2008 R2 (64-bit) | Microsoft Corporation | | | |
| Name | Publisher | | | | | | | | | | |
| Coffalyser - Microsoft SQL Server 2008 R2 (Preconfig... | MRC-Holland | | | | | | | | | | |
| Coffalyser.Net | MRC-Holland | | | | | | | | | | |
| Microsoft SQL Server 2008 R2 (64-bit) | Microsoft Corporation | | | | | | | | | | |
| 5.  Click Yes to confirm you want to uninstall Coffalyser.Net | | | | | | | | | | | |





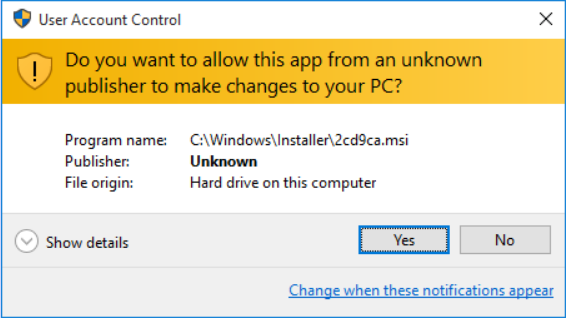
| | | |
|--|---|---|
| 6. |  | |
| If prompted, click Yes in the User Account Control window | | |
| 7. | Close the Control Panel of Windows |  |
| 8. | Proceed with the installation of the new version of Coffalyser.Net as described in Procedure installing new version – preconfigured SQL server or Procedure installing new version – not preconfigured SQL server | |

PROCEDURE INSTALLING NEW VERSION – PRECONFIGURED SQL SERVER

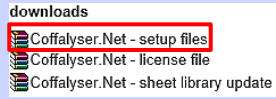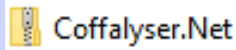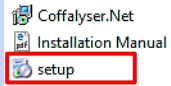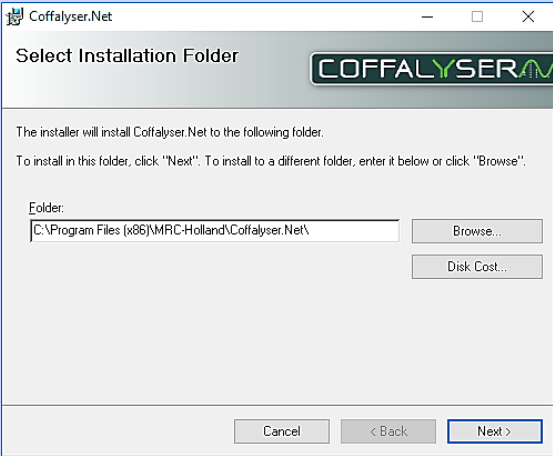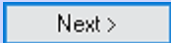
This procedure should be followed when for the original installation an SQL database server was used that was preconfigured for Coffalyser.Net. In case for the original installation a previously installed, not preconfigured SQL database server was used, see **Procedure installing new version – not preconfigured SQL server**

| | |
|--|---|
| <p>1. Download the zip file with the Coffalyser.Net – setup files from your MyMLPA account on www.mlpa.com</p> |  |
| <p>2. Extract the zip file</p> |  |
| <p>3. Start the installation by opening the setup file.</p> |  |
| <p>The Coffalyser.Net installation wizard opens</p> | |
| <p>4. Maximize the Coffalyser.Net installation wizard from the task bar if this doesn't happen automatically.</p> | |
| <p>5.</p>  <p>If it is desired to install Coffalyser.Net on another location than the default, click Browse. If the default location is ok, continue with the next step</p> | |
| <p>6. Click Next</p> |  |
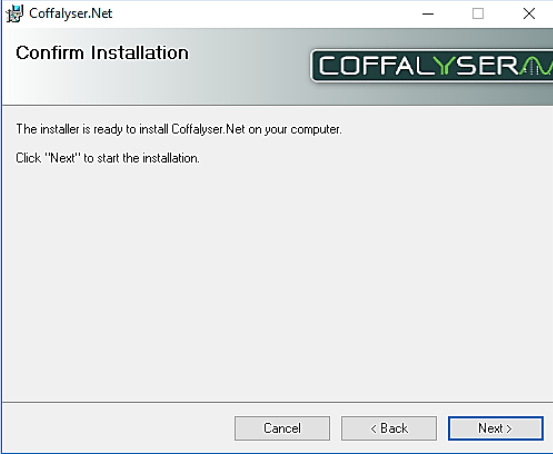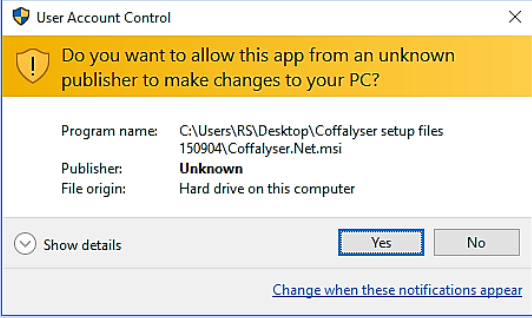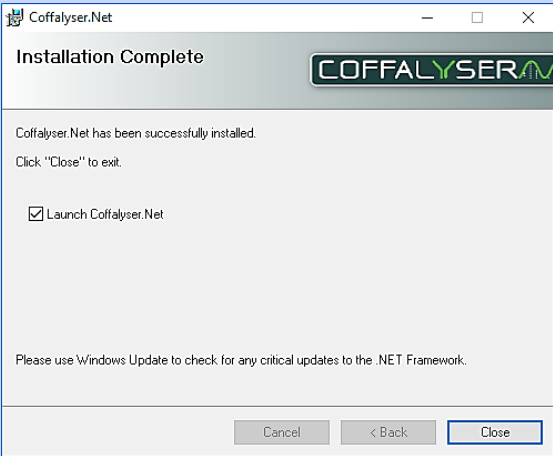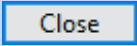
| | | |
|-----|---|---|
| 7. |  | |
| 8. | <p>Click Next to start the installation</p>  | |
| 9. | <p>If prompted, click Yes in the User Account Control window</p>  | |
| 10. | <p>Ensure that the selection box <i>Launch Coffalyser.Net</i> is checked</p> <p>Click Close to exit the installation window</p> |  |
| | <p>Coffalyser.Net starts</p> | |

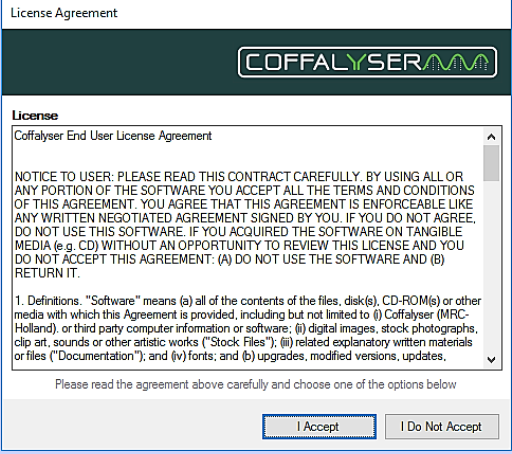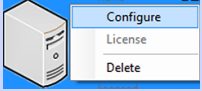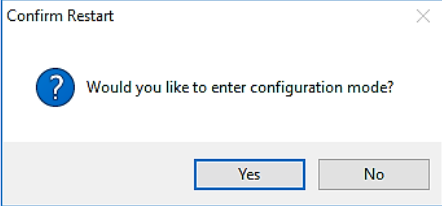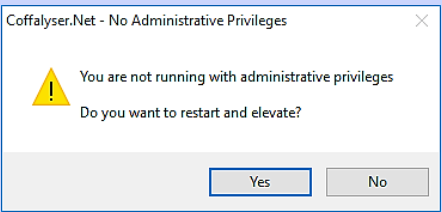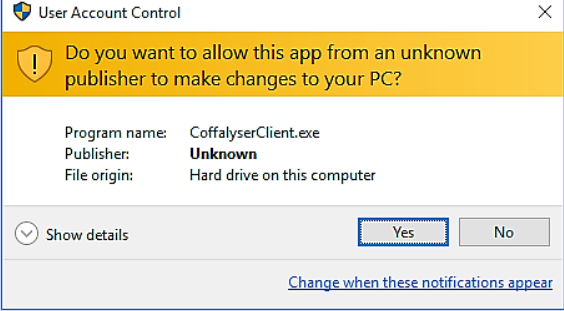
| | | |
|--|--|---|
| 11. |  <p>License Agreement</p> <p>Coffalyser</p> <p>License</p> <p>Coffalyser End User License Agreement</p> <p>NOTICE TO USER: PLEASE READ THIS CONTRACT CAREFULLY. BY USING ALL OR ANY PORTION OF THE SOFTWARE YOU ACCEPT ALL THE TERMS AND CONDITIONS OF THIS AGREEMENT. YOU AGREE THAT THIS AGREEMENT IS ENFORCEABLE LIKE ANY WRITTEN NEGOTIATED AGREEMENT SIGNED BY YOU. IF YOU DO NOT AGREE, DO NOT USE THIS SOFTWARE. IF YOU ACQUIRED THE SOFTWARE ON TANGIBLE MEDIA (e.g. CD) WITHOUT AN OPPORTUNITY TO REVIEW THIS LICENSE AND YOU DO NOT ACCEPT THIS AGREEMENT: (A) DO NOT USE THE SOFTWARE AND (B) RETURN IT.</p> <p>1. Definitions: "Software" means (a) all of the contents of the files, disk(s), CD-ROM(s) or other media with which this Agreement is provided, including but not limited to (i) Coffalyser (MRC-Holland), or third party computer information or software; (ii) digital images, stock photographs, clip art, sounds or other artistic works ("Stock Files"); (iii) related explanatory written materials or files ("Documentation"); and (iv) fonts; and (b) upgrades, modified versions, updates.</p> <p>Please read the agreement above carefully and choose one of the options below</p> <p><input type="button" value="I Accept"/> <input type="button" value="I Do Not Accept"/></p> | |
| 12. | <p>In the Coffalyser.Net Server Selection window right mouse click on the (grey) server with the name of your computer and select <i>Configure</i></p> |  |
| 13. |  <p>Confirm Restart</p> <p>Would you like to enter configuration mode?</p> <p><input type="button" value="Yes"/> <input type="button" value="No"/></p> | |
| 14. |  <p>Coffalyser.Net - No Administrative Privileges</p> <p>You are not running with administrative privileges Do you want to restart and elevate?</p> <p><input type="button" value="Yes"/> <input type="button" value="No"/></p> | |
| 15. |  <p>User Account Control</p> <p>Do you want to allow this app from an unknown publisher to make changes to your PC?</p> <p>Program name: CoffalyserClient.exe Publisher: Unknown File origin: Hard drive on this computer</p> <p><input type="button" value="Yes"/> <input type="button" value="No"/></p> <p>Show details</p> <p>Change when these notifications appear</p> | |
| <p>If prompted, click Yes in the User Account Control window</p> <p>The Coffalyser.Net – Configuration Wizard opens</p> | | |

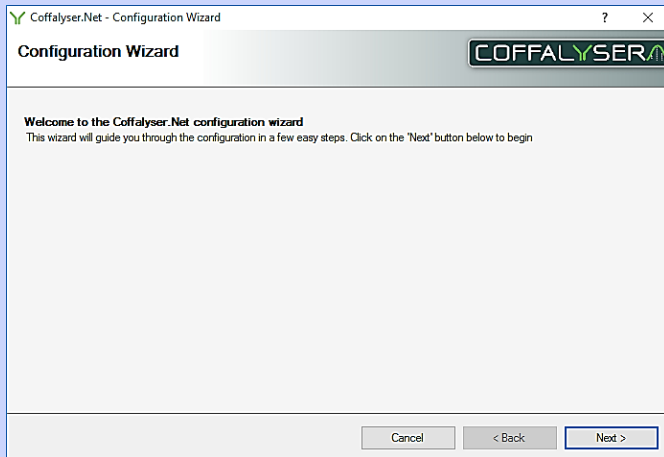16.



Click **Next**

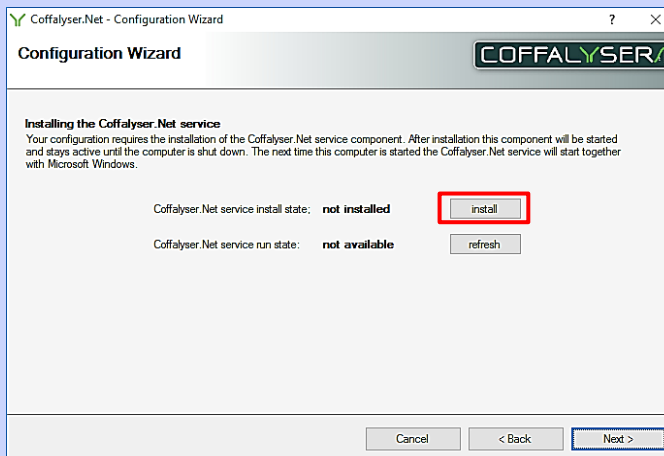17.




Click **install** to install the Coffalyser.Net service

18.




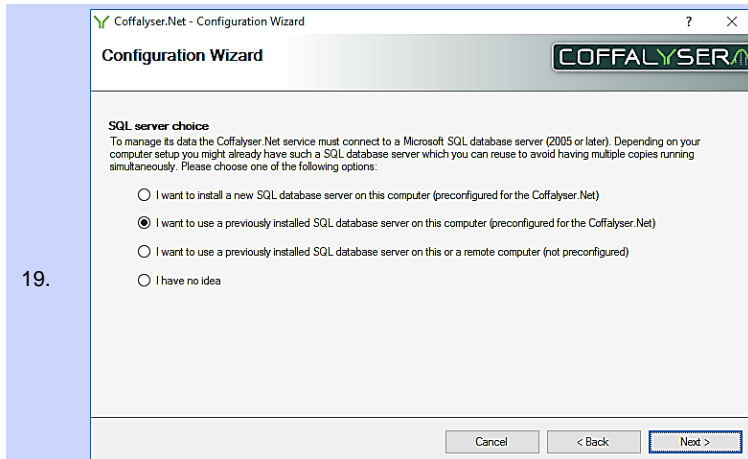
Click **Next**



19.

Select the option *I want to use a previously installed SQL database server on this computer (preconfigured for the Coffalyser.Net)* if it is not preselected

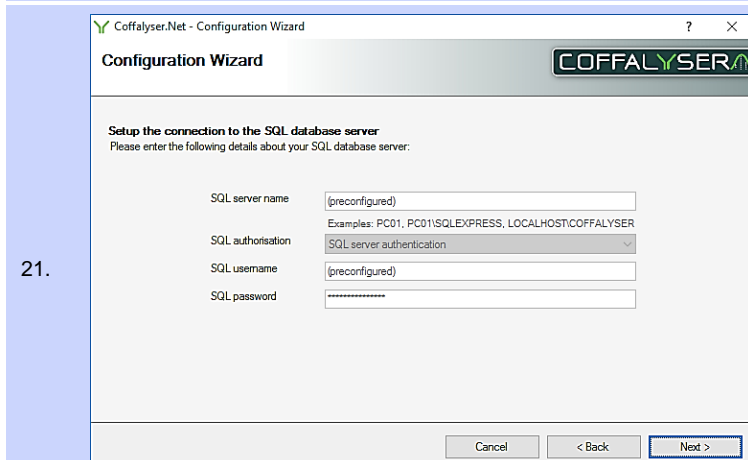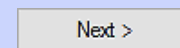20. Click **Next**



21.

Click **Next**

Do not change the database name

22.

Coffalyser.Net - Configuration Wizard

Configuration Wizard COFFALYSER

Database name
Please enter the name of the Coffalyser.Net database

database name

Cancel < Back Next >

Click **Next**

23.

Coffalyser.Net - Configuration Wizard

Configuration Wizard COFFALYSER

Upgrading the existing Coffalyser.Net database
The database you selected is outdated. It must be upgraded before it can be used in combination with the installed Coffalyser.Net service. Be aware that upgrading the database might out date other Coffalyser installations, which might have to be upgraded as well.

upgrade the database

Cancel < Back Next >

If prompted, check the selection box *upgrade the database*

24. Click **Next**

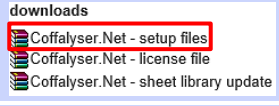
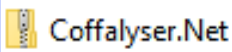
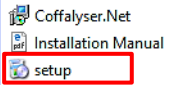
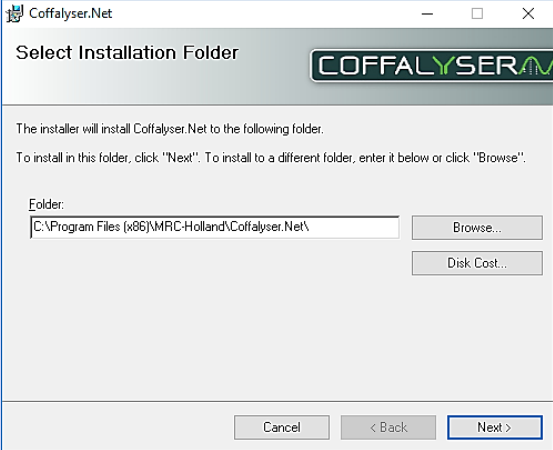
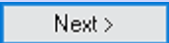
Next >

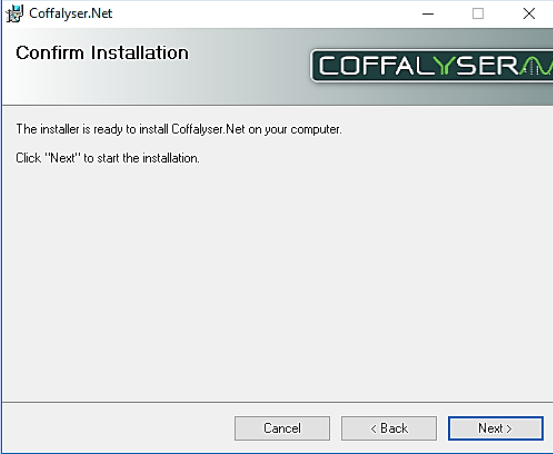
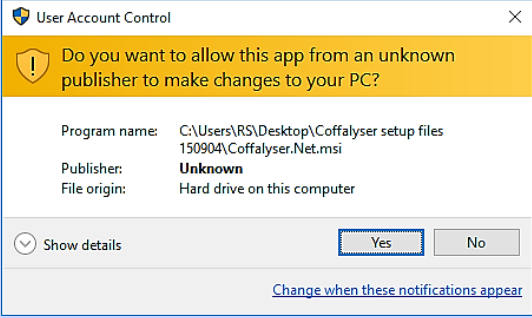
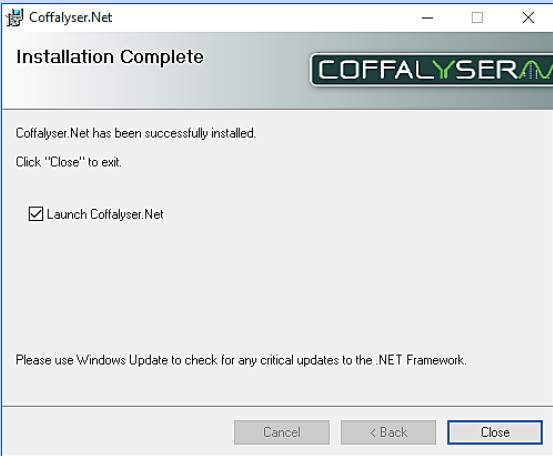
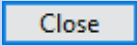
Do not change the server description


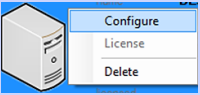
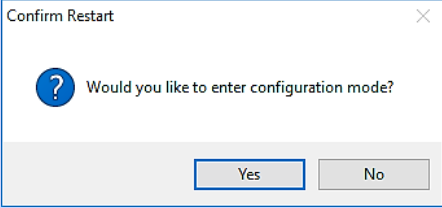
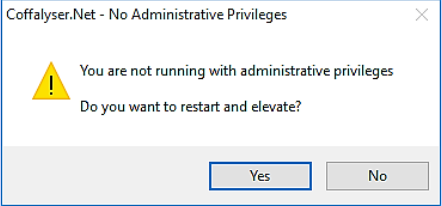
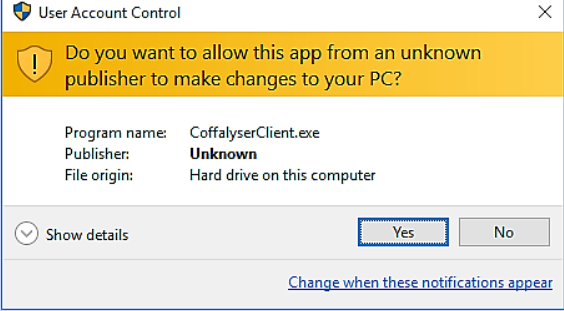
| | | |
|-----|---|---|
| 25. |  <p>Click Finish</p> | |
| 26. |  <p>Ensure that the selection box <i>start the Coffalyser after the wizard is closed</i> is checked</p> | |
| 27. | Click Close to exit the Configuration wizard |  |
| 28. | If the server in the Coffalyser.Net Server Selection window is orange, install your license as described in the chapter License on page 23 | |
| 29. | Coffalyser.Net is now ready for use | |

PROCEDURE INSTALLING NEW VERSION – NOT PRECONFIGURED SQL SERVER


This procedure should be followed when for the original installation a previously installed, not preconfigured SQL database server was used. In case for the original installation an SQL database server was used that was preconfigured for Coffalyser.Net., see **Procedure installing new version – preconfigured SQL server**

| | |
|---|---|
| 1. Download the zip file with the Coffalyser.Net – setup files from your MyMLPA account on www.mlpa.com |  |
| 2. Extract the zip file |  |
| 3. Start the installation by opening the setup file. |  |
| The Coffalyser.Net installation wizard opens | |
| 4.  <p>If it is desired to install Coffalyser.Net on another location than the default, click Browse. If the default location is ok, continue with the next step</p> | |
| 5. Click Next |  |

| | | |
|-----------------------|--|---|
| 6. |  | |
| 7. | <p>Click Next to start the installation</p>  | |
| 8. | <p>If prompted, click Yes in the User Account Control window</p>  <p>Ensure that the selection box <i>Launch Coffalyser.Net</i> is checked</p> | |
| 9. | <p>Click Close to exit the installation window</p> |  |
| Coffalyser.Net starts | | |


| | | |
|--|--|---|
| 10. |  <p>Click I Accept to accept the license agreement of Coffalyser.Net</p> | |
| 11. | <p>In the Coffalyser.Net Server Selection window <u>right</u> mouse click on the (grey) server with the name of your computer and select <i>Configure</i></p> |  |
| 12. |  <p>Click Yes to enter the configuration mode</p> | |
| 13. |  <p>If prompted, click Yes to restart and elevate in order to run with administrative privileges</p> | |
| 14. |  <p>If prompted, click Yes in the User Account Control window</p> | |
| <p>The Coffalyser.Net – Configuration Wizard opens</p> | | |

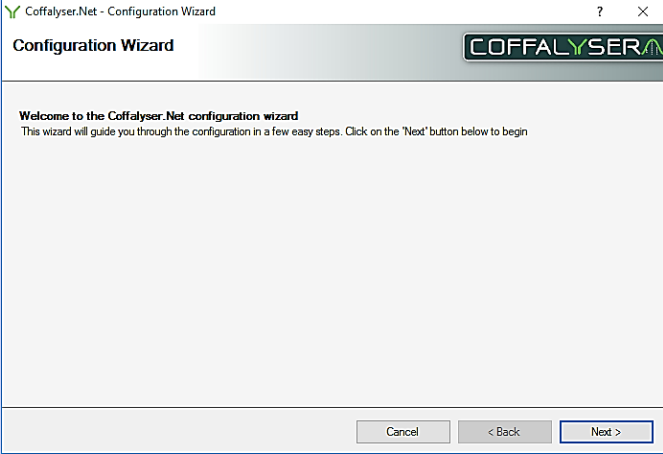15.




Click **Next**

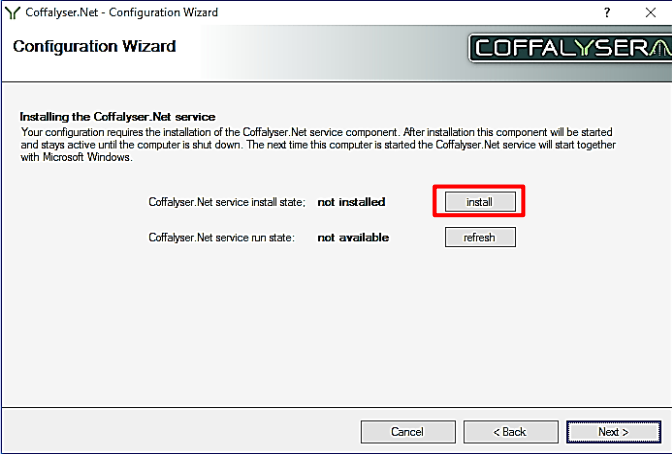16.




Click **install** to install the Coffalyser.Net service

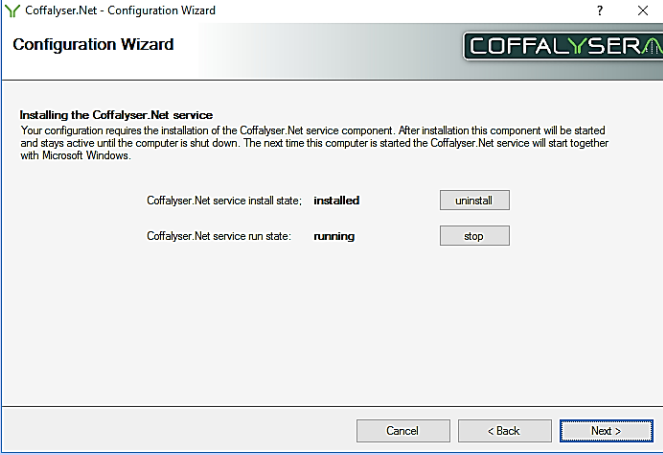17.





Click **Next**

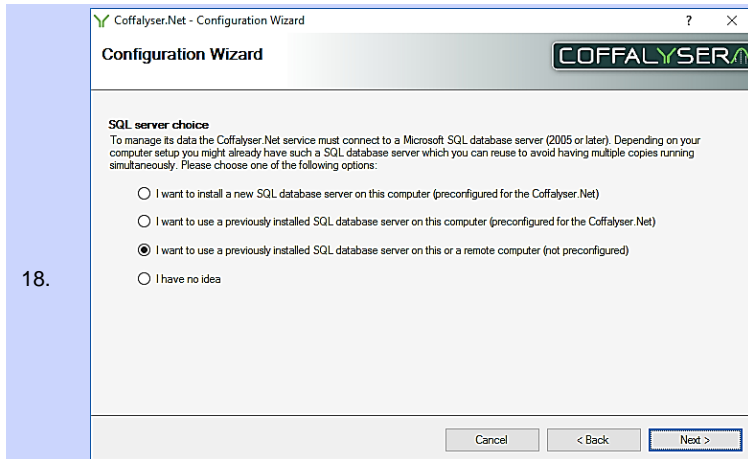

18.

Select the option *I want to use a previously installed SQL database server on this or a remote computer (not preconfigured)* if it is not preselected

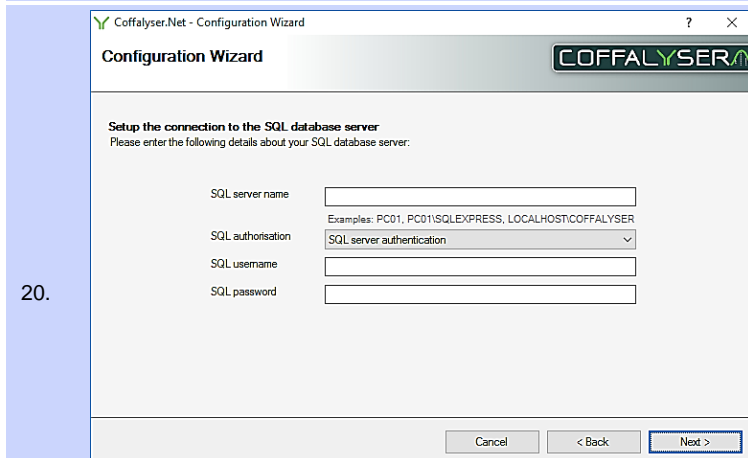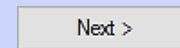
19. Click **Next**




20.

Enter the information of the SQL server in the designated fields if it is not filled in yet

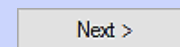
21. Click **Next**



Do not change the database name

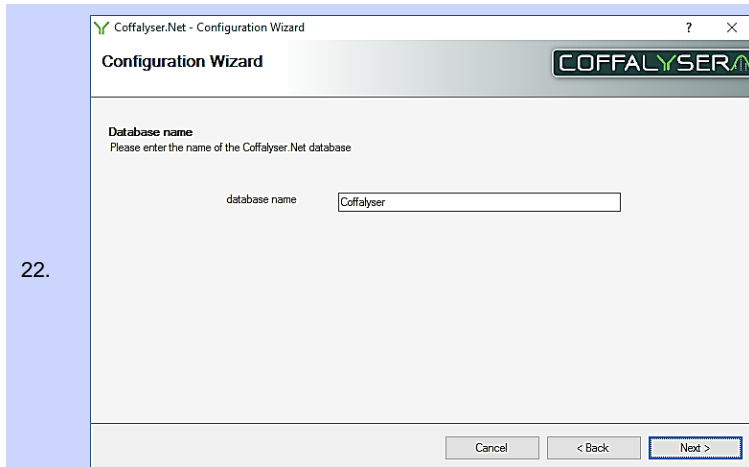


Click **Next**

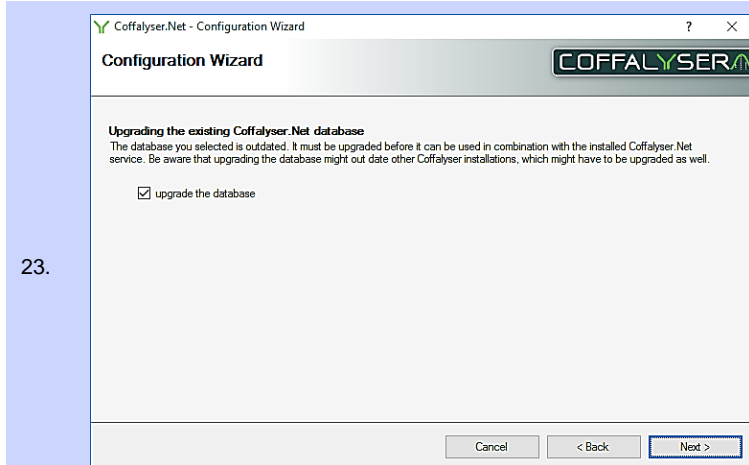

If prompted, check the selection box *upgrade the database*

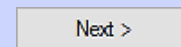24. Click **Next**




Do not change the server description

| | | |
|-----|---|---|
| 25. |  <p>Click Finish</p> | |
| 26. |  <p>Ensure that the selection box <i>start the Coffalyser after the wizard is closed</i> is checked</p> | |
| 27. | Click Close to exit the Configuration wizard |  |
| 28. | If the server in the Coffalyser.Net Server Selection window is orange, install your license as described in the chapter License on page 23 | |
| 29. | Coffalyser.Net is now ready for use | |

More information & contact

More information about Coffalyser.Net can be found in the Coffalyser.Net Reference Manual which is available online at www.mlpa.com.

| | |
|---|---|
|  | |
| | MRC-Holland BV; Willem Schoutenstraat 1 1057 DL, Amsterdam, The Netherlands |
| Website | www.mlpa.com ; www.mlpa.eu |
| E-mail | info@mlpa.com (information & technical questions); order@mlpa.com (orders) |
| Phone | +31 888 657 200 |