

# MRC Holland Support

Support > Help Centre > MLPA & Coffalyser.Net > Coffalyser.Net Data Analysis Software > Data Analysis > Comparative Analysis > [How can I export all sample reports at once in Coffalyser.Net?](#)

---


## How can I export all sample reports at once in Coffalyser.Net?

This article was retrieved from [support.mrcholland.com](https://support.mrcholland.com) on Friday, 9th May 2025.

### Note

This article is part of the [Coffalyser.Net Tips & Tricks](#) collection.

Coffalyser.Net conveniently allows you to export sample reports of all samples simultaneously.

- In the comparative analysis tab , right-click on a sample and select Open Sample Results.
- You can now start the export of all sample reports in three ways:
  1. Navigate to the Sample result overview or Sample report tab, right-click and select Create PDF Reports All Samples.
  2. In the Sample result overview or Sample report tab, press the keyboard shortcut Ctrl + Alt + R.
  3. In any tab, click the  button in the toolbar.
    - In the dialog box that opens, enter a name for the folder in which the reports will be saved in the File name field, and select the location where you want this folder to be created.
- Click Save to create the folder and sample reports.

Comparative Analysis Sample Results Explorer

127-REF-P081-B1-1109-KTU-DNA2 [R]

140-REF-P081-B1-1109-KTU-DNA2 [R]

161-REF-P081-B1-1109-KTU-DNA2 [R]

183-REF-P081-B1-1109-KTU-DNA2 [R]

R0145-sample-P081-B1-1109-KTU-DNA

Sample result overview

Sample report

ratio chart

Electropherograms

127-REF-P081-B1-1109-K

Height

Area

NF1-1

17-026.445712

2920

29010

NF1-2

17-026.445712

2920

29010

NF1-4

17-026.445712

2920

29010

NF1-6

17-026.445712

2920

29010

NF1-11

17-026.445712

2920

29010

NF1-13

17-026.445712

2920

29010

NF1-15

17-026.445712

2920

29010

NF1-17

17-026.445712

2920

29010

NF1-18

17-026.445712

2920

29010

NF1-21

17-026.445712

2920

29010

NF1-23

17-026.445712

2920

29010

NF1-26

17-026.445712

2920

29010

NF1-28

17-026.445712

2920

29010

NF1-29

17-026.445712

2920

29010

NF1-32

17-026.445712

2920

29010

NF1-35

17-026.445712

2920

29010

OMG-2

17-026.445712

2920

29010

NF1-37

17-026.445712

2920

29010

NF1-39

17-026.445712

2920

29010

NF1-41

17-026.687519

3533

39825

17q (n=27)

Export Grid Data

Ctrl+E

3300

Export Grid Data All Samples

Ctrl+Alt+E

4512

Export Grid To Image

Ctrl+S

4048

Export Grid To Image All Samples

Ctrl+Alt+S

8384

Create PDF Report

Ctrl+R

5100

Create PDF Reports All Samples

Ctrl+Alt+R

2237

Region Analysis

0757

Create separate reference region

8088

Resize Column Width

0337

Resize Row Height

9860

Sort Data

2842

Hide Column

9600

Show All Columns

5008

Channels

0298

Copy Data To Clipboard

9962

Copy To Clipboard

0782

Tags

Coffalyser.Net

## Disclaimer

The information provided in this material is correct for the majority of our products. However, for certain applications, the instructions for use may differ. In the event of conflicting information, the relevant instructions for use take precedence.