

MRC Holland Support

Support > Help Centre > MLPA & Coffalyser.Net > Coffalyser.Net Data Analysis Software > Data Analysis > Comparative Analysis > [How can I export all sample reports at once in Coffalyser.Net?](#)


How can I export all sample reports at once in Coffalyser.Net?

This article was retrieved from support.mrcholland.com on Saturday, 27th April 2024.

Note

This article is part of the [Coffalyser.Net Tips & Tricks](#) collection.

Coffalyser.Net conveniently allows you to export sample reports of all samples simultaneously.

- In the comparative analysis tab , right-click on a sample and select Open Sample Results.
- You can now start the export of all sample reports in three ways:
 1. Navigate to the Sample result overview or Sample report tab, right-click and select Create PDF Reports All Samples.
 2. In the Sample result overview or Sample report tab, press the keyboard shortcut Ctrl + Alt + R.
 3. In any tab, click the  button in the toolbar.
 - In the dialog box that opens, enter a name for the folder in which the reports will be saved in the File name field, and select the location where you want this folder to be created.
- Click Save to create the folder and sample reports.

