

# MRC Holland Support

Support > Help Centre > MLPA & Coffalyser.Net > Coffalyser.Net Data Analysis Software > Data Analysis > Comparative Analysis > How can I export all sample reports at once in Coffalyser.Net?

---


## How can I export all sample reports at once in Coffalyser.Net?

This article was retrieved from [support.mrcholland.com](https://support.mrcholland.com) on Monday, 2nd October 2023.

### Note

This article is part of the [Coffalyser.Net Tips & Tricks](#) collection.

Coffalyser.Net conveniently allows you to export sample reports of all samples simultaneously.

- In the comparative analysis tab , right-click on a sample and select Open Sample Results.
- You can now start the export of all sample reports in three ways:
  1. Navigate to the Sample result overview or Sample report tab, right-click and select Create PDF Reports All Samples.
  2. In the Sample result overview or Sample report tab, press the keyboard shortcut Ctrl + Alt + R.
  3. In any tab, click the  button in the toolbar.
    - In the dialog box that opens, enter a name for the folder in which the reports will be saved in the File name field, and select the location where you want this folder to be created.
- Click Save to create the folder and sample reports.

Comparative Analysis Sample Results Explorer

Sample result overview | Sample report | ratio chart | Electropherograms

127-REF-P081-B1-1109-KTU-DNA2 [R]  
 140-REF-P081-B1-1109-KTU-DNA2 [R]  
 161-REF-P081-B1-1109-KTU-DNA2 [R]  
 183-REF-P081-B1-1109-KTU-DNA2 [R]  
 R0145-sample-P081-B1-1109-KTU-DNA

Probe target info		Height	Area
NF1-1	17-026.445712	2920	29010
NF1-2	17-026.445712		0300
NF1-4	17-026.445712		4512
NF1-6	17-026.445712		4048
NF1-11	17-026.445712		8384
NF1-13	17-026.445712		5100
NF1-15	17-026.445712		2237
NF1-17	17-026.445712		5620
NF1-18	17-026.445712		0757
NF1-21	17-026.445712		8088
NF1-23	17-026.445712		0337
NF1-26	17-026.445712		9860
NF1-28	17-026.445712		2842
NF1-29	17-026.445712		9600
NF1-32	17-026.445712		5008
NF1-35	17-026.445712		0298
OMG-2	17-026.445712		3846
NF1-37	17-026.445712		9962
NF1-39	17-026.445712		0782
NF1-41	17-026.687519	3533	39825
NF1-42	17-026.687519	1331	17025

17q (n=27)

- Export Grid Data Ctrl+E
- Export Grid Data All Samples Ctrl+Alt+E
- Export Grid To Image Ctrl+S
- Export Grid To Image All Samples Ctrl+Alt+S
- Create PDF Report Ctrl+R
- Create PDF Reports All Samples Ctrl+Alt+R**
- Region Analysis
- Create separate reference region
- Resize Column Width
- Resize Row Height
- Sort Data
- Hide Column
- Show All Columns
- Channels
- Copy Data To Clipboard
- Copy To Clipboard

Tags  
Coffalyser.Net

### Disclaimer

*The information provided in this material is correct for the majority of our products. However, for certain applications, the instructions for use may differ. In the event of conflicting information, the relevant instructions for use take precedence.*