

MRC Holland Support

Support > Help Centre > MLPA & Coffalyser.Net > Coffalyser.Net Data Analysis Software > Tips & Tricks for Coffalyser.Net > [How can I resize data grids in Coffalyser.Net?](#)

How can I resize data grids in Coffalyser.Net?



This article was retrieved from support.mrcholland.com on Sunday, 11th May 2025.

Note



This article is part of the [Coffalyser.Net Tips & Tricks](#) collection.

You can resize the rows and columns of all data grids in order to fit them on your screen. This is particularly useful for experiments with a large number of probes and/or a large number of samples.

Column widths can be adjusted in the following ways:

- With the   buttons in the toolbar above the grids.
- With the keyboard shortcuts Ctrl + Shift + Minus and Ctrl + Shift + Plus.
- By clicking the right mouse button and selecting the desired option from the Resize Column Width submenu.

Similarly, row heights can be adjusted in the following ways:

- With the   buttons in the toolbar above the grids.
- With the keyboard shortcuts Ctrl + Alt + Minus and Ctrl + Alt + Plus.
- By clicking the right mouse button and selecting the desired option from the Resize Row Height submenu.

Tags

Coffalyser.Net

Related Pages

- [How can I zoom in and out in Coffalyser.Net?](#)

Disclaimer

The information provided in this material is correct for the majority of our products. However, for certain applications, the instructions for use may differ. In the event of conflicting information, the relevant instructions for use take precedence.