

MRC Holland Support

Support > Help Centre > digitalMLPA & Coffalyser digitalMLPA > Coffalyser digitalMLPA Data Analysis Software > Installation & Configuration > [How can I update Coffalyser digitalMLPA to the latest available version?](#)

How can I update Coffalyser digitalMLPA to the latest available version?

This article was retrieved from support.mrcholland.com on Sunday, 11th May 2025.

Note

This article is about Coffalyser digitalMLPA for digitalMLPA data analysis. For information about Coffalyser.Net for conventional MLPA data analysis, see [this article](#). Updating Coffalyser digitalMLPA is a very straightforward procedure.

1. Remove the old folder with Coffalyser digitalMLPA from your system, or move it to an archive location.
2. Log in to your account on our website.
3. [Download](#) the latest Coffalyser digitalMLPA software package.
4. Extract the latest Coffalyser digitalMLPA software package to a folder on your hard drive.

Security policies or antivirus software may prevent running downloaded software from a network location. If this happens, we recommend moving the package to a local directory (e.g. *C:\Coffalyser digitalMLPA*).

Tags

Coffalyser digitalMLPA

Related Files

- [Coffalyser digitalMLPA User Manual](#)

Disclaimer

The information provided in this material is correct for the majority of our products. However, for certain applications, the instructions for use may differ. In the event of conflicting information, the relevant instructions for use take precedence.