

# MRC Holland Support

Support > Help Centre > MLPA & Coffalyser.Net > Coffalyser.Net Data Analysis Software > Data Analysis > Sheet Library & CE Devices > [How do I add a CE device to Coffalyser.Net?](#)

---

## How do I add a CE device to Coffalyser.Net?

This article was retrieved from [support.mrcholland.com](https://support.mrcholland.com) on Sunday, 11th May 2025.

A capillary electrophoresis (CE) device has to be added to your organisation before you can analyse MLPA data with Coffalyser.Net. You can add a CE device as follows:

1. Right-click on CE Devices and select Add CE Device ....
2. On the general tab, select the CE device from the CE device dropdown menu.
3. Select the filter set from the CE device filter dropdown menu.
4. Optionally enter something in the location and remarks fields to recognize the CE device.
5. Click OK.

This procedure is described in more detail in Appendix V of the [reference manual](#). Alternatively, you can watch the video below.

In case you are unable to view the video on YouTube you can also watch it [here](#).

## Background

Each type of capillary electrophoresis device has its own specifications, which is why a specific CE device has to be selected for each experiment. Appendix V of the [reference manual](#) contains more information on the supported devices and filter sets. The appendix also contains instructions on how to view, edit and delete a CE device in Coffalyser.Net.

Tags

Coffalyser.Net

Instruction Videos

---

### Related Pages

- [Video: How to analyse SALSA MS-MLPA data in Coffalyser.Net](#)
- [Video: How to analyse SALSA MLPA data in Coffalyser.Net](#)

---

Disclaimer

*The information provided in this material is correct for the majority of our products. However, for certain applications, the instructions for use may differ. In the event of conflicting information, the relevant instructions for use take precedence.*