

# MRC Holland Support

Support > Help Centre > Coffalyser.Net > Installation & Configuration > Licenses > How do I find the email address associated with my Coffalyser.Net license?


## How do I find the email address associated with my Coffalyser.Net license?

This article was retrieved from support.mrcholland.com on Tuesday, 9th August 2022.

Every Coffalyser.Net server requires a free license associated with an email address. It is important to ensure that this email address is current, as we need to be able to contact you in case of important (safety) notifications about the software. More information about our licenses can be found [here](#).

If you are not sure which email address was used for your license, you can easily find out as follows:

1. Start Coffalyser.Net.
2. In the *Coffalyser.Net Server Selection* dialog, right-click on the server and select *License*. If Coffalyser.Net starts with the *Login* dialog instead, you can bring up the server selection dialog by clicking on the *select* button next to the server address.

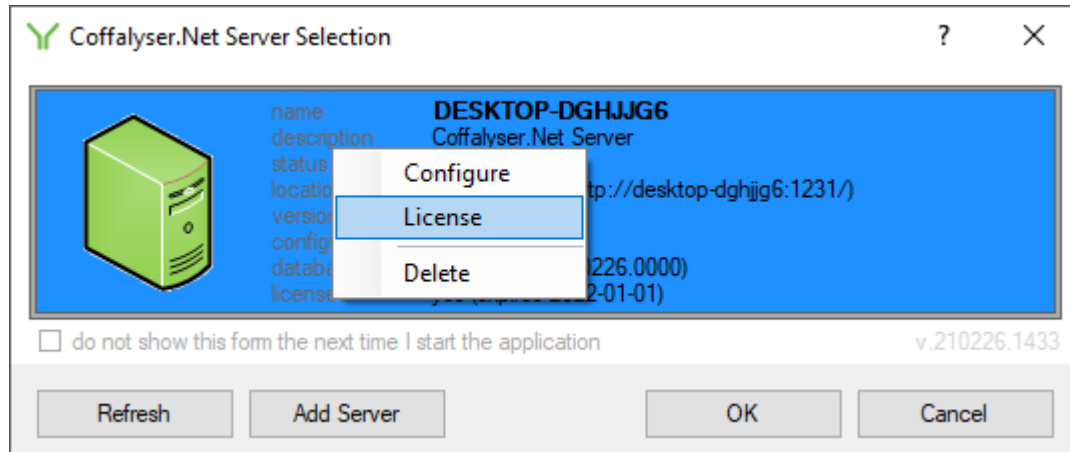

3. The dialog that opens shows your license, the email address associated with the license, and when the license expires ([read how to renew your license](#)).

You can also:

1. Start Coffalyser.Net.
2. Log in as you normally would.
3. Click on *About* in the menu. Your license information will be listed in the *License* section.

[Contact us](#) if you need help changing your email address, or if you no longer have access to the email address associated with your license.

Tags

Coffalyser.Net

#### Related Pages

---

- [Why do Coffalyser.Net and Coffalyser digitalMLPA require a license?](#)

#### Disclaimer

---

*The information provided in this material is correct for the majority of our products. However, for certain applications, the instructions for use may differ. In the event of conflicting information, the relevant instructions for use take precedence.*