

MRC Holland Support

Support > Help Centre > Coffalyser.Net > Working with Coffalyser.Net > Fragment Analysis > How do I interpret the probe counter in Coffalyser.Net?




How do I interpret the probe counter in Coffalyser.Net?

This article was retrieved from support.mrcholland.com on Tuesday, 9th August 2022.

The probe counter in Coffalyser.Net provides a visual warning when an unexpected number of probes have been found in a sample. The probe counter has the following format:

number of probes found / number of probes expected | number of probes found that were *not* expected

The last component is only shown if there were unexpected probe signals.



probes	
	50/50
	49/50
	50/50 1

It is important to inspect the data whenever the probe counter provides a warning, as an unexpected count can be caused by biological properties of the sample but also by experimental or data analysis issues.

Background

Probes for which a signal is expected



In many probemixes, all probes are expected to have a signal in healthy individuals. If a lower number of probes is found, this may be caused by a homozygous deletion, but also by e.g. an issue with peak identification ([binning](#)) or signals that are too low. The example below shows a probemix with 50 probes that are expected to produce a signal. Signals were found for all 50 probes in the first sample, but only for 49 probes in the second sample.

probes	
	50/50
	49/50

Probes for which a signal is *not* expected











Some probemixes have probes for which a signal is *not* expected in healthy individuals, such as mutation-specific probes. If a significant signal for a mutation-specific probe is found, this is unexpected and will result in a probe counter such as for the second sample in the screenshot below. The data should be inspected to

determine if the sample might be positive for the mutation or if the result was caused by an artefact such as a nonspecific peak.













probes
 50/50
 50/50 1

Differences between samples in the number of probes expected

The number of probes expected may differ between samples in the same experiment. If the probemix contains one or more probes for the Y chromosome, a signal for these probes is expected in male samples but not in female samples. In the example below there is one probe for the Y chromosome.

probes	DNA	DD	X	Y
 57/57				
 56/56				

In MS-MLPA, the number of expected probes in digested and undigested samples may differ because some probes may be expected to produce no signal after digestion in healthy individuals. If the probemix has one or more digestion control probes, these are also expected to be absent after digestion. The example below is for a probemix in which eight probes are not expected to give a signal in the digested reaction. In the last digested sample, six of these nevertheless gave a signal. This may be due to the methylation status of the sample DNA, but could also be caused by e.g. incomplete digestion or nonspecific peaks.

sample type	bin smpl	FRSS	FMRS	probes
sample	<input type="checkbox"/>			 27/27
digested sample	<input type="checkbox"/>			 19/19
sample	<input type="checkbox"/>			 27/27
digested sample	<input type="checkbox"/>			 19/19 6

Criteria for a probe to be counted in the probe counter

Probes are only counted as 'found' when their signal is:

- Above the minimum peak amplitude for probes (in RFU) defined for the electrophoresis device.
- Above 10% of the median signal of the reference probes.

Probes that are below 10% of the median signal of the reference probes but above the device's minimum signal may still be identified and present in the results, but will not be counted in the probe counter. This helps to more easily identify unexpected signals using the probe counter.

In rare cases, slope correction may influence whether a probe signal crosses the

abovementioned thresholds. As a result, the probe counter may change after the comparative analysis, when slope correction is performed. If this happens there were probably one or more probes with a signal just around the thresholds for inclusion in the probe counter.

Warnings for unexpected probe counts

An orange warning is displayed for an unexpected probe count. Test samples also get a subtraction to their FMRS score if a *reference probe* is missing, as this is very unusual. If reference samples are missing the signal of any expected probe, a red warning is displayed for the probe counter and the FMRS score gets a heavy penalty, because samples that miss an expected probe signal are not suitable as reference samples. In the example below, the second *reference* sample and the last *test* sample both miss a signal for a target probe, but this results in different warnings.

sample type	bin smpl	FRSS	FMRS	probes
reference	<input type="checkbox"/>			50/50
reference	<input type="checkbox"/>			49/50
sample	<input type="checkbox"/>			50/50
sample	<input type="checkbox"/>			49/50

The information in this article is for Coffalyser.Net [v.210226.1433](#) and newer. In Coffalyser.Net v.140721.1958, the probe counter did not yet include the clear distinction between expected and unexpected probe signals found.

Tags

Coffalyser.Net

Related Pages

- [What do the probe colours in the electropherogram and genomic profile in Coffalyser.Net mean?](#)

Disclaimer

The information provided in this material is correct for the majority of our products. However, for certain applications, the instructions for use may differ. In the event of conflicting information, the relevant instructions for use take precedence.