

MRC Holland Support

Support > Help Centre > MLPA & Coffalyser.Net > Coffalyser.Net Data Analysis Software > Error Messages > [How do I fix SQL Server errors while installing Coffalyser.Net?](#)

How do I fix SQL Server errors while installing Coffalyser.Net?

This article was retrieved from support.mrcholland.com on Saturday, 20th April 2024.

Installation and setup of Microsoft SQL Server may fail if one of the [system requirements](#) are not met. Check if the requirements are met, and retry the installation if possible.

In some environments, installation of the preconfigured SQL Server may be problematic for a variety of reasons. In such cases, advanced users are recommended to manually install a newer version of Microsoft SQL Server, or to use an existing SQL Server installation if possible. Coffalyser.Net works with all editions of SQL Server that are currently supported by Microsoft, and can connect to SQL Servers that are hosted elsewhere in the network if needed.

Please [contact us](#) for personal assistance if you keep experiencing issues while installing SQL Server – in our experience, remaining issues are usually very specific to a particular system, and personalized support is often required.

Background

Coffalyser.Net uses a Microsoft SQL Server database to store all data. As part of the installation procedure of Coffalyser.Net, Microsoft SQL Server 2022 is downloaded and installed. Coffalyser.Net usually installs without issues, but in some cases there are compatibility issues between Microsoft SQL Server and specific settings or properties of a computer.

Tags

Coffalyser.Net

Related Pages

- [What are the system requirements for Coffalyser.Net?](#)

Disclaimer

The information provided in this material is correct for the majority of our products.

However, for certain applications, the instructions for use may differ. In the event of conflicting information, the relevant instructions for use take precedence.